**AGIMATIC-S  
AGIMATIC-N  
AGIMATIC-E**

**7000270  
7000271  
7000273**

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## GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) ***IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.***
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.*

## PACKING LIST

The standard equipment consist of the following components:

DESCRIPTION	CODE	QUANTITY
AGIMATIC S	7000270	1
AGIMATIC N	7000271	1
AGIMATIC E	7000273	1
Power supply wire	7140	1
Safety ring	2049	1
Teflon magnet 8Øx39mm	1000039	1
Instruction manual	80002	1

## ACCESSORIES

<b>Denomination</b>	<b>Code</b>
<i>Magnetic stirring-bar 7Øx9mm</i>	1000009
<i>Magnetic stirring-bar 4Øx12mm</i>	1000012
<i>Magnetic stirring-bar 7Øx19mm</i>	1000019
<i>Magnetic stirring-bar 8Øx32mm</i>	1000032
<i>Magnetic stirring-bar 8Øx39mm</i>	1000039
<i>Magnetic stirring-bar 10Øx45mm</i>	1000045
<i>Magnetic stirring-bar 12Øx75mm</i>	1001061
<i>Magnetic stirring-bar 16Øx102mm</i>	1001062
<i>Magnetic stirring-bar 16Øx127mm</i>	1001063
<i>Magnetic stirring-bar 19Øx150mm</i>	1001064
<i>Triangular magnetic stirring-bar TRIMAG 10 (sidex35 length)</i>	1000536
<i>Triangular magnetic stirring-bar TRIMAG 14 (sidex39 length)</i>	1000537
<i>Triangular magnetic stirring-bar TRIMAG 14 (sidex45 length)</i>	1000538
<i>Triangular magnetic stirring-bar TRIMAG 14 (sidex55 length)</i>	1000539
<i>Magnetic stirring-bar ROTOR-PAT 10Øx32 mm</i>	1000001
<i>OVAL magnetic stirring-bar 16 Øx30 mm length</i>	1000788
<i>OVAL magnetic stirring-bar 20 Øx17 mm length</i>	1000789
<i>Crosshead magnetic stirring bars 25Øx15 mm height</i>	1000790

<i>Floating magnet IMANFLOTER. Structure 50 mm. Magnet 39 mm</i>	1000897
<i>Floating magnet IMANFLOTER. Structure 60 mm. Magnet 45 mm</i>	1000898
<i>Concentric magnet CENTRIMAN (length 32 mm)</i>	1000900
<i>Concentric magnet CENTRIMAN (length 50 mm)</i>	1000901
<i>Concentric magnet CENTRIMAN (length 50 mm)</i>	1000901
<i>Concentric magnet CENTRIMAN (length 55 mm)</i>	1000902
<i>Stirring bar retriever</i>	1000020
<i>Box with assorted magnetic stirring bars</i>	1000487
<i>Pt100 probe with handle which allows read and control liquid temperature (model E)</i>	1001610
<i>Digital contact thermometer SENSOTERM -40-270°C (models N &amp; E)</i>	1001009
<i>Electric contact thermometer 0-100°C (models N &amp; E)</i>	3521159
<i>Electric contact thermometer 0-200°C (models N &amp; E)</i>	3522159
<i>Electric contact thermometer 0-300°C (models N &amp; E)</i>	3523159
<i>Stay rod 12Øx500mm length</i>	1000725
<i>Clamp to hold thermometer</i>	7000025
<i>Stainless steel round tray 2 litres</i>	1001059
<i>Reducing discs set with cover</i>	3000917
<i>Bosshead</i>	1000069

## **EQUIPMENT DESCRIPTION**

*The magnetic electronic stirrers described in this manual comprise three models with stainless steel plate.*

*The stirrers **AGIMATIC S, N and E** is a group of stirrers with analogic technology for the control of speed and temperature. This technology allows setting parameters in a precise and repetible way.*

*S model allows digital reading and control of speed.*

*N model allows setting a heating level as a rate of total heating power. To regulate the temperature of the stirred liquid in a better way, use an electronic controller Sensoterm II as an accessory.*

*E model allows setting the heater plate temperature. Including 1001610 probe as an accessory, it allows controlling liquid temperature at same same time that it is stirring.*

*In some application it is important bearing in mind that heater plate temperature and the liquid one are very different.*

POWER	AGIMATIC S	25W
	AGIMATIC N	550W
	AGIMATIC E	
Maximum temperature (N & E models)		350°C
Temperatura resolution (E model)		± 2°C
Heating control (N & E models)		N:0...100% E:50...350°C
External temperature control connection (N & E models)		Yes
Speed		60...1600rpm
Stirring power		10.000ml
Dimensions		14,5x19x26cm
Weight (S model)		3,0Kg
Weight (N & E models)		3,2Kg
Temperatura limiter (N & E models)		370...390°C

## INSTALLATION

*Place the stirrer on a flat, horizontal, level surface near the mains.*

*Connect the connection cable at the rear part of the stirrer base and then plug it into the mains.*

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*Be sure that the voltage supply is the same as the one indicated on the characteristics plate.*




*Do not use the equipment if it is not earthed.*

*If you change the plug, bear in mind the following:*

<b>Blue cable:</b>	<b>Neutral</b>
<b>Brown cable:</b>	<b>Phase</b>
<b>Yellow/green cable:</b>	<b>Earth</b>

## OPERATION

### CONTROL PANEL: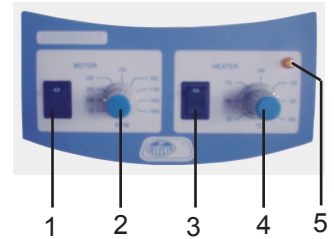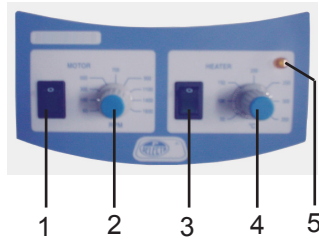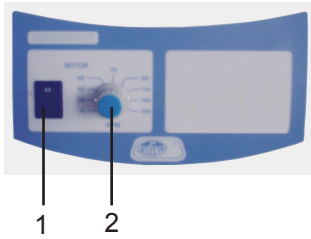

#### S MODEL




1. Agitation switch
2. Speed set button

#### N MODEL

1. Agitation switch
2. Speed set button
3. Heating switch
4. Set % heating power
5. Heating indicator lamp

#### E MODEL

1. Agitation switch
2. Speed set button
3. Heating switch
4. Set temperature
5. Heating indicator lamp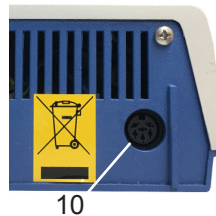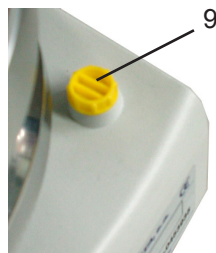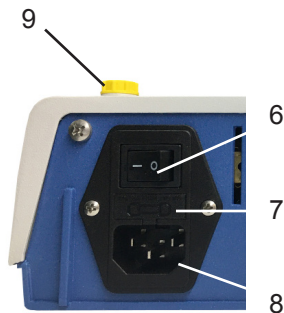

#### REAR PART

6. Main switch
7. Fuse carrier
8. Mains plug
9. Thermometer support rod housing
10. External control connection or liquid probe (N & E models)

**STARTING UP :****AGIMATIC-S:**

Once the indications in the **Installation** section have been made, switch on the main switch on the rear part, place the recipient to be stirred on the top plate, let the teflon-coated magnetic stirring bar drop into the recipient, turn on the "MOTOR" switch and adjust speed with "RPM" knob.

**AGIMATIC-N:**

For its starting up operation and to select the speed, do as the above model, AGIMATIC-S. There is no heating until motor is working.

Then switch on the HEATER switch, and with the regulator for heating power, select the required "%". The yellow lamp will show the working of the heater.

**AGIMATIC-E:**

To start it working and to adjust speed, follow the above indications. With the regulator "HEATER" the suitable working temperature may be selected with the "°C" knob.

**LIQUID TEMPERATURE CONTROL**

(N & E models)

1. Connection to work without external connection of a temperature control device.

2. Connection for contact thermometer.

- Put the thermometer with the bulb submerged into the liquid. Connect the thermometer in the stirrer rear part plug.
- Set the temperature turning the superior part of the thermometer. The stirrer must be configured to work at maximum of power rate or temperature.
- The connection is as shown on figure 1.

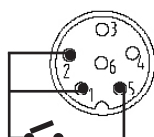
Figure 1

3. Connection for regulator "SENSOTERM II" (1001009)

- Put the probe submerged into the liquid. Connect the probe in the stirrer rear part plug.
- Set the temperature on the «SENSOTERM II».
- Set the power rate or temperature on the stirrer at maximum.
- The connection is as shown on figure 2.

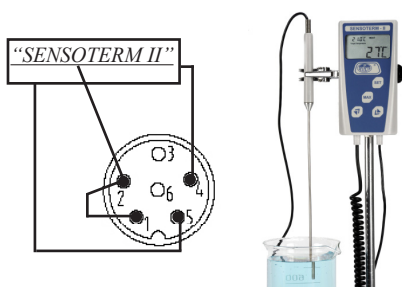
Figure 2

## MAINTENANCE

**Before removing the stirrer casing, disconnect the equipment from the mains.**

**The guarantee will not have any effect if the equipment is manipulated by somebody who has not been expressly authorized by J.P.SELECTA, s.a.u.**

### CLEANING:

For the cleaning of the different parts, use the following products:

Cleaning of stainless steel: Alcohol

Cleaning of plastic: Alcohol with a cotton duster.

## WARRANTY

*This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA,s.a.u.*

*Any manipulation of the equipment by not authorized personnel by J.P. SELECTA,s.a.u. automatically cancels the guarantee.*

## SPARE PARTS

Heating plate	0032026
Control circuit:	
for model S	0038515
for model N	0038514
for model E	0038513
Pattern magnet	0015449
Motor	0024283
Pt100 probe	0043028 (model E only)