PRECISTERM 2L.	6000137 / 6000385
PRECISTERM 5L.	6000138 / 6000386
PRECISTERM 12L.	6000140 / 6000387
PRECISTERM 20L.	6000141 / 6000388
PRECISTERM 45L.	6000142 / 6000389

BAÑOS TERMOSTÁTICOS PARA AGUA Y ACEITE
THERMOSTATIC BATHS FOR WATER AND OIL

INFORMACIÓN GENERAL

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

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PACKING LIST

The standard equipment consist of the following components:

Descripción / Description	Código / Code	Cant. / Qty
Precistern 2L.	6000137	1
Precistern 2L. 200°C	6000385	1
Precistern 5L.	6000138	1
Precistern 5L. 200°C	6000386	1
Precistern 12L.	6000140	1
Precistern 12L. 200°C	6000387	1
Precistern 20L.	6000141	1
Precistern 20L. 200°C	6000388	1
Precistern 45L.	6000142	1
Precistern 45L. 200°C	6000389	1
Manual de instrucciones / Instruction manual	80003	1

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ACCESSORIES

(The following components, at user's request, must be fitted in factory)

Denominación	Código	Denomination	Code
NIVEL DE QUITA Y PON Para baños que deban usarse sin tapa de disco.	6001400	REMOVABLE LEVEL For baths to be used without perforated cover.	6001400
NIVEL FIJO Debe acoplarse en fábrica.	6000123	FIXED LEVEL It must be fitted in factory.	6000123
TERMOMETRO LECTOR Para acoplar con nuez (120°C) Para acoplar con nuez (200°C)	1000120 1000200	THERMOMETER To adapt with bosshead (120°C) To adapt with bosshead (200°C)	1000120 1000200
TERMOMETRO LECTOR Para acoplar a tapa tejadillo (120°C) Para acoplar a tapa tejadillo (200°C)	1000872 1000873	THERMOMETER To be fitted at gable cover (120°C) To be fitted at gable cover (200°C)	1000872 1000873
NUEZ SOPORTE PARA TERMÓMETRO LECTOR Abertura de fijación hasta 40mm. En acero inoxidable 304.	6000896	SUPPORT BOSSHEAD FOR THERMOMETER Set opening up to 40mm . In AISI 304 stainless steel.	6000896

Para baño Código For bath code		Tapa con orificios y discos reductores de 105, 80, 60 y 37mm Ø, con alojamiento para termómetro lector. Cover with holes and reducing discs 105, 80, 60 and 37 mm dia., with a hole for the thermometer in AISI 304 stainless steel	Tapa Tejadillo Gable cover in AISI 304 stainless steel	Bandeja cubrerresistencias Perforated shelf in AISI 304 stainless steel
6000137	6000385	6003231 (1 hole)	No	6000011
6000138	6000386	6003232 (2 holes)	6000710	6000022
6000140	6000387	6003234 (4 holes)	6000711	6000044
6000141	6000388	6003236 (6 holes)	6000712	6000066
6000142*	6000389*	2x6003236 (6 holes) = 12 holes)	6000714	6000088

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*Because of the large surface of baths code 6000142 and 6000389, to work at temperatures above 50°C, the use of cover with holes and reducing discs or gable cover is recommended

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Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Código Code	Temperatura máxima °C Maximum temperature °C	Capacidad litros Capacity litres	Medidas útiles cm. Tank dimensions cm			Medidas exteriores cm. Overall dimensions cm			Consumo W Consumption W	Peso Kg Weight Kg
			Alto Depth	Ancho Length	Fondo Width	Alto Depth	Ancho Length	Fondo Width		
6000137	110	2	15	15	13	20	20	27	390	3.5
6000385	200								640	
6000138	110	5	15	30	14	20	37	27	600	6
6000386	200								1110	
6000140	110	12	15	30	30	20	40	42	1110	8
6000387	200								1400	
6000141	110	20	15	48	30	20	58	42	1500	10
6000388	200								2100	
6000142*	110	45	15	62	50	22	69	62	2400	18
6000389*	200								3800	

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The temperature stability is ±1 K for baths of 100°C and ±2 K for baths of 200°C.

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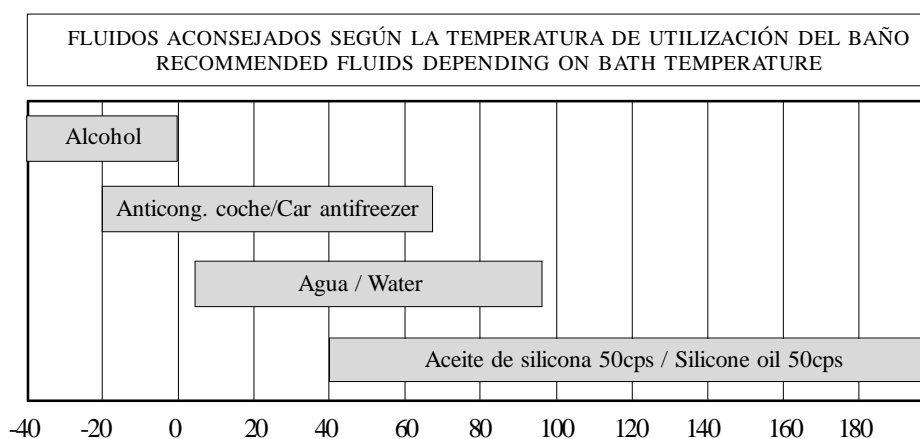
EQUIPMENT DESCRIPTION

PRECISTERM baths with adjustable liquid temperature from ambient +5°C up to 110°C (Stability of ±1°C) or up to 200°C (Stability of ±2°C), depending on the model. To use the baths fill the tank with liquid until the level covers at least the heater element. The liquid or mixture to be employed must be adequated to the model of bath and the working temperature. As a guidance, here are some liquids (to use these liquids, follow the manufacturer's instructions):

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INSTALLATION

Place the bath on a flat, horizontal and level surface. Bear in mind the weight of the liquid when choosing the support. Provide for bath drainage (bath 6000142 and 6000389).

Be sure that the voltage supply is the same as the one indicated on the characteristics plate.

Do not use the apparatus if it is not earthed.

If you change the plug bear in mind the following:

***Blue cable: Neutral.
Brown cable: Phase.
Yellow/green cable: Earth.***

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It is recommended the use of decalcified water. DO NOT USE DEIONIZED OR DEMINERALIZED WATER as it has a great tendency to rust. To prolong the life of the bath and avoid rusting of the parts, dissolve 1 gr of CO₃Na₂ (Sodium Carbonate) per litre of water used in the bath.

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
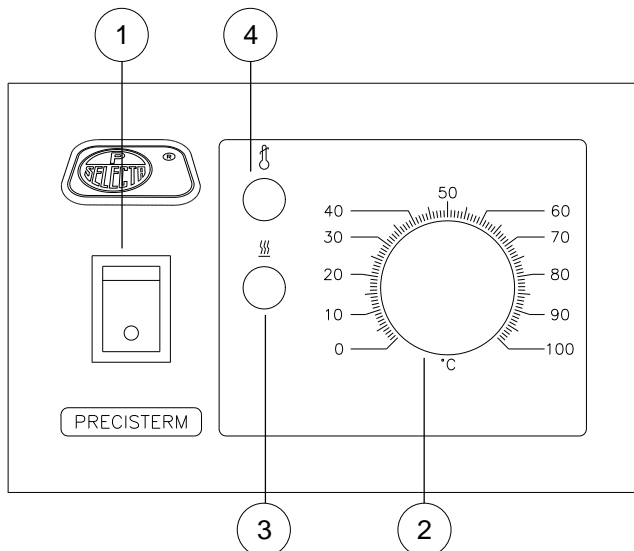
OPERATION

CONTROL PANEL:

1. Main switch with indicator lamp.
2. Temperature control knob.
3. Heater operation indicator lamp (amber).
4. Safety thermostat operation indicator lamp (red).

REAR PART

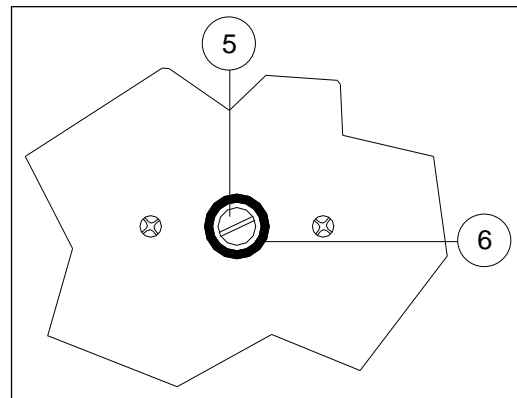5. Safety thermostat control screw.
6. Safety thermostat manual reset button (grey).

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STARTING UP OPERATION :

1. Fill the bath with liquid. Leave at least 3 cm without filling to avoid splashes.
2. Connect the bath to the mains.
3. Turn on the main switch (1). The Green indicator lamp will light up.
4. Turn on the temperature control thermostat (2) from low to high according to the working temperature. The heater operation indicator lamp (amber) will light up (3).
5. After a certain time, the heater operation indicator lamp (3) will begin to operate intermittently. When the cycle of these intermittences becomes uniform, that will indicate that temperature of liquid is stabilized.

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6. *The thermostat scale graduation is only useful as a reference, because the real temperature is the one shown by a thermometer introduced into the liquid.*
7. *Be careful with liquid evaporation. The heater element must ALWAYS be covered with liquid. Replenish when necessary.*

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ADJUSTMENT FOR THE SAFETY THERMOSTAT:

The low limit of working temperature for the safety thermostat is nearly 50°C. The desired temperature of the safety thermostat must be at least 5°C above the working temperature. The adjusting process of the safety thermostat is long. For that reason, the safety thermostat is useful when the users work habitually at the same temperature.

1. *Turn the safety thermostat screw (5) clockwise to maximum.*
2. *Select with the temperature control knob (2) the desired temperature of the safety thermostat.*
3. *Turn on the bath and leave it to stabilize.*
4. *When the bath is stabilized, slowly turn the safety thermostat screw (5) anticlockwise until it clicks and the red indicator lamp (4) lights up.*
5. *Wait until the bath cools to reset the safety thermostat with (6) button.*
6. *Select the working temperature of the bath with temperature control knob (2).*

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SPARE PARTS


To guarantee the safety of the equipment, the spare parts must be purchased from J.P.SELECTA, S.A.

DESCRIPCIÓN / DESCRIPTION	CÓDIGO / CODE									
Modelo / Model	6000137	6000385	6000138	6000386	6000140	6000387	6000141	6000388	6000142	6000389
1. Resistencia Heating element	39063	39049	39061	39050	39051	39060	39012	39059	39123	39124
2. Termostato de seguridad Safety thermostat	43016	43017	43016	43017	43016	43017	43016	43017	43016	43017
3. Termostato de regulación Control thermostat	43006	43005	43006	43005	43006	43005	43006	43005	43006	43005
4. Racord termostato + tuerca termostato Thermostat racord + thermostat nut	34141 + 34142									
5. Cable manguera / Mains cable	7024									
6. Contactador Omron / Omron contactor	No									13012

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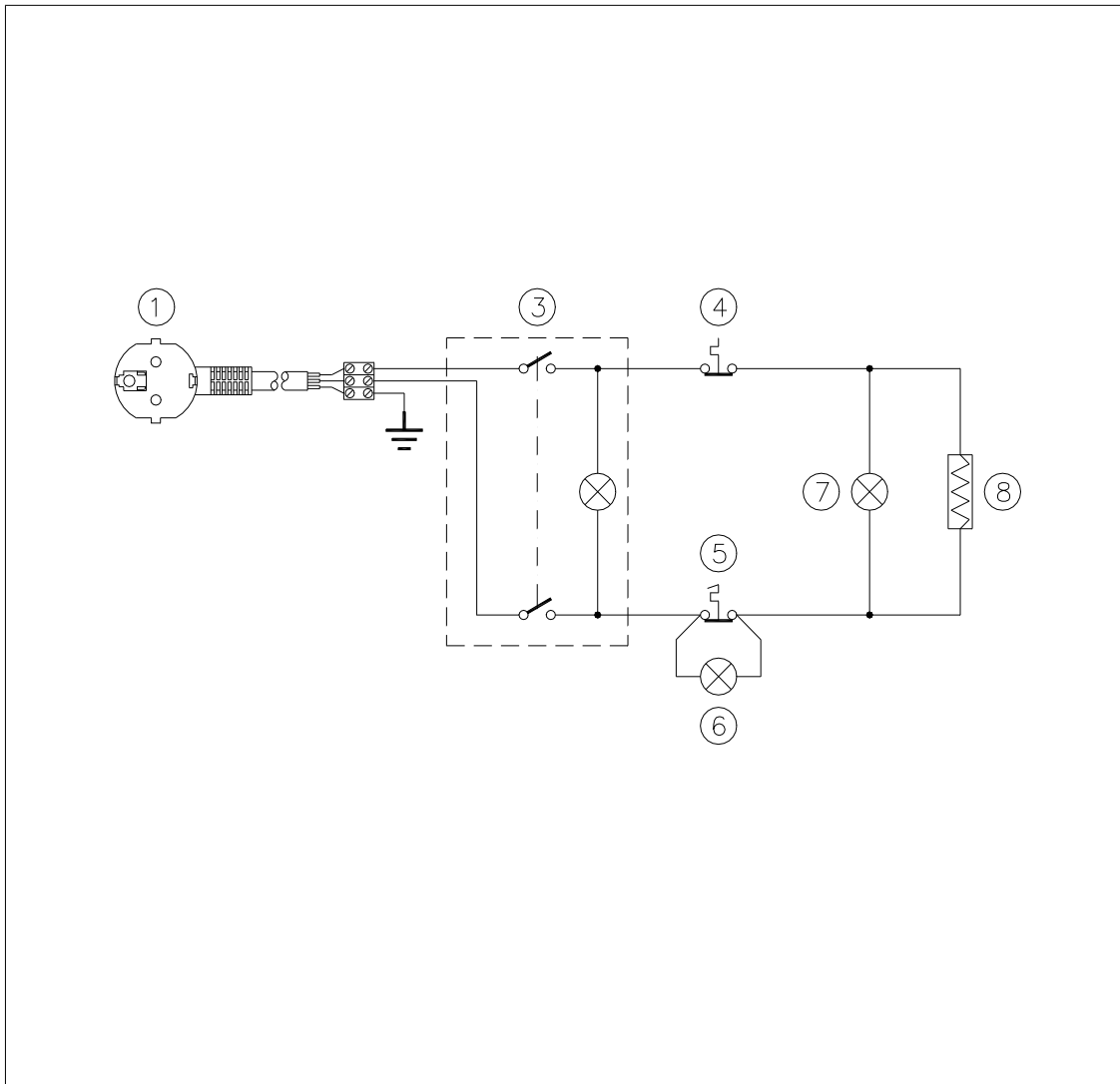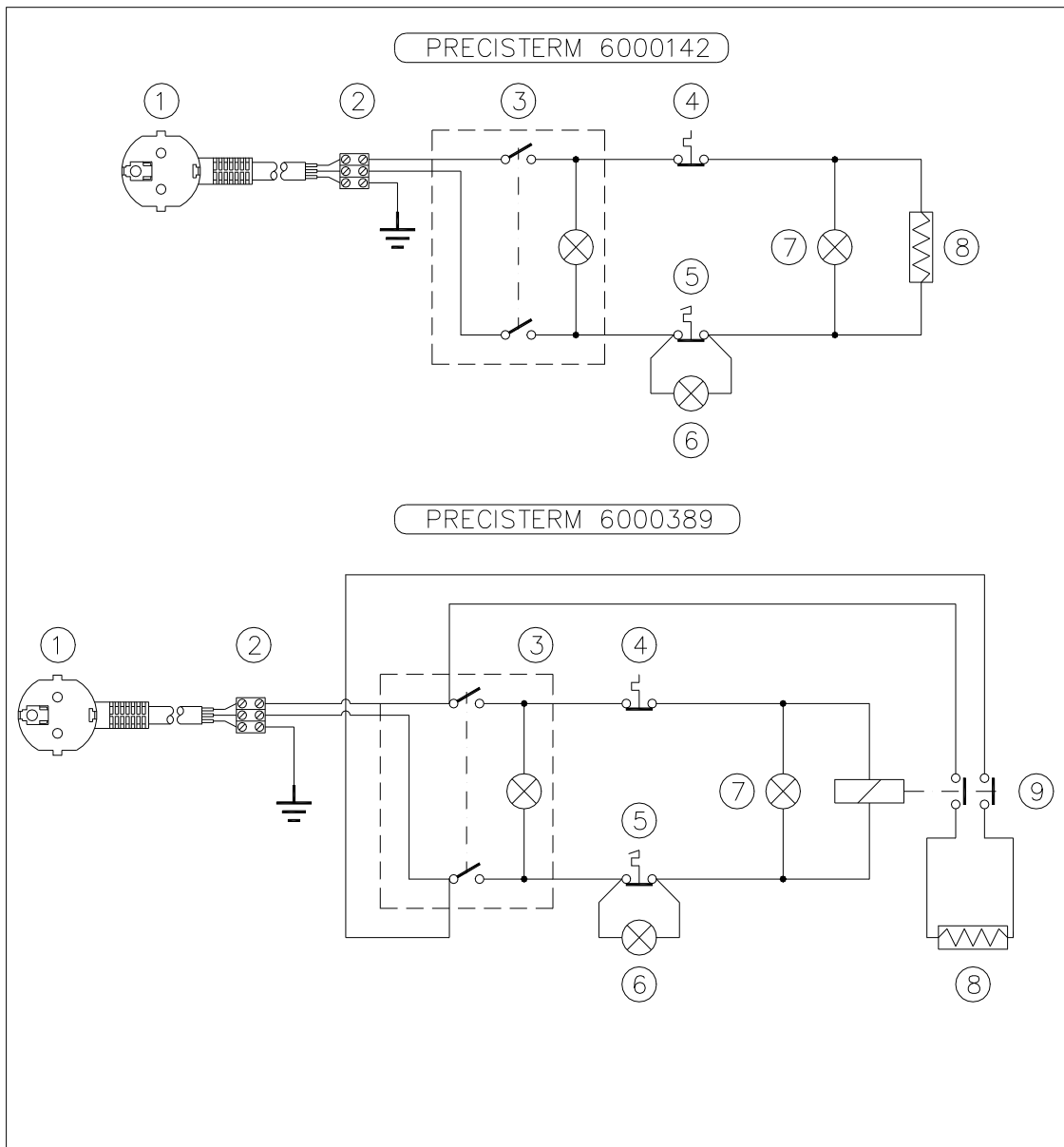
MAINTENANCE

Before removing the cover disconnect the apparatus from the mains.

The control elements and heating elements must only be manipulated by authorized personnel.

The distilled or demineralized water has a great tendency to rust. To prolong the life of the bath and avoid rusting of the parts, dissolve 1 gr of CO_3Na_2 (Carbonate of Soda) per litre of water used in the bath.

CLEANING:

If normal tapwater is used calcareous deposits will appear in the bath and in the heating element. A heater covered with lime should not be used as it can lose heating power and deteriorate more quickly. Periodical cleaning of the inside of the tank and the heating element with a solution of 10% of 28% Chlorhydric acid or a solution of water and vinegar (Acetic acid) is advised. Rinse with clean water afterwards.

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: alcohol

Cleaning of plastic: Alcohol with cotton duster.

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use or causes beyond the control of J.P. SELECTA, S.A.

Any manipulation of the apparatus by personnel not authorized by J.P. SELECTA, S.A. automatically cancels the guarantee.

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