AGIMAN

7000272

**AGITADOR MAGNÉTICO PARA MANTAS CALEFACTORAS
MAGNETIC STIRRER FOR HEATING MANTLES**

INDICE

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) ***IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.***
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.*

PACKING LIST

The standard equipment consists of the following components:

Piece	Code	Quantity
<i>Magnetic stirrer AGIMAN</i>	7000272	1
<i>Attachment device for heating mantles</i>		1
<i>Rotor-Pat magnetic stirring bar 10Øx32</i>	1000001	1
<i>Instruction manual</i>	80008	1

ACCESSORIES

<i>Denomination</i>	<i>Code</i>
<i>PTFE-COATED STIRRING BARS</i>	
<i>Magnetic stirring bar Rotor-Pat 10Ø máx.x32mm length.</i>	1000001
<i>Oval magnetic stirring bar 16Øx30mm length.</i>	1000788
<i>Oval magnetic stirring bar 20Øx17mm length.</i>	1000789
<i>STIRRING BAR RETRIEVER</i>	1000020
<i>Flexible, PTFE-coated. Length: 280mm</i>	
<i>CONTACT THERMOMETER</i>	
<i>Range: 0-100°C Scale division: 1°C</i>	3521159
<i>Range: 0-200°C Scale division: 2°C</i>	3522159
<i>Range: 0-300°C Scale division: 2°C</i>	3523159
<i>Total length: 395mm. Immersion stem length: 150mm</i>	
<i>SENSOTERM II digital contact thermometer</i>	
<i>Range -50°C to 300°C Resolution: 1°C</i>	1001009
<i>STAY ROD</i>	1000725
<i>Ø12mm x 500mm length.</i>	
<i>CLAMP TO HOLD THERMOMETER</i>	7000025
<i>BOSSHEAD</i>	1000069
<i>HEATING MANTLES</i>	3000447
<i>Flask capacity 250ml. Consumption 130W</i>	
<i>HEATING MANTLES</i>	3000448
<i>Flask capacity 500ml. Consumption 270W</i>	
<i>HEATING MANTLES</i>	3000449
<i>Flask capacity 1000ml. Consumption 410W</i>	
<i>HEATING MANTLES</i>	3000450
<i>Flask capacity 2000ml. Consumption 530W</i>	

TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

<i>Code</i>	<i>Overall dimensions cm.</i>			<i>Consumption W.</i>	<i>Weight Kg.</i>
	<i>Height</i>	<i>Length</i>	<i>Width</i>		
7000272	14,5	19	26	25	3,5

EQUIPMENT DESCRIPTION

Magnetic stirrers **AGIMAN** are designed for adjustable temperatures up to 300°C with an electronic regulator of heat power synchronised with heater-on indicator lamp. Also the agitation speed can be adjusted up to 1600 r.p.m. with an electronic speed regulator.

INSTALLATION

During the first starting up, you will notice that the heating mantle smokes and has a characteristic smell. It is due to the newness of the heating material. These circumstances are completely normal and will disappear after a short time.

For a correct operation and a better performance, it is recommended the use of adjustable flasks that match the mantle capacity.

Avoid spilling liquids on the heating material surface, as it would immediately destroy it.

CAUTION!!! IMPORTANT FOR YOUR SAFETY

Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the equipment.


Do not use the equipment if it is NOT earthed.

If you have to change the plug, bear in mind the following:

Blue cable:	Neutral.
Brown cable:	Phase.
Yellow/green cable:	Earth.

OPERATION

CONTROL PANEL:

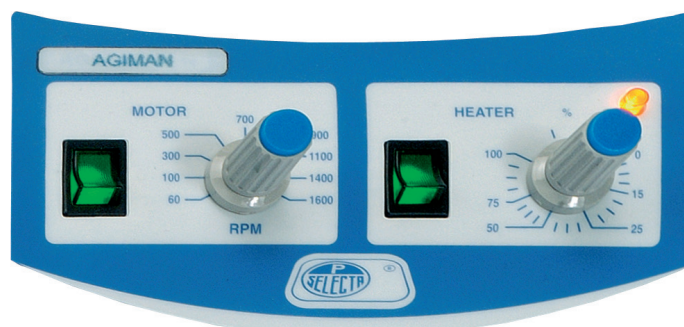To place the heating mantle in the stirrer, a circular perforated plate must be used and its plug must be connected to the base of the socket at the back of the stirrer.

Once all the preceding indications are checked, place the recipient to be stirred on the stirrer plate, let the PTFE-coated magnetic stirring bar drop into the recipient, turn on the "MOTOR" switch and adjust the speed with "RPM" knob.

For working with heating, switch on the "HEATER" switch, and select the required "%" with the heating power regulator.

TEMPERATURE REGULATION WITH CONTACT THERMOMETER (optional accessory):

Place the thermometer with the bulb submerged into the liquid, supported by a special clamp (code 7000025), a bosshead (code 1000069) and a stay rod (code 1000725). Connect the thermometer plug at the rear part of the equipment. Adjust the working temperature by turning the thermometer regulator and fixing it conveniently.

Once the temperature has been selected, turn on the "HEATER" switch and place the selector "%" at the maximum position. Start up the stirrer, adjust the working speed and turn on the "HEATER" switch.

MAINTENANCE

Before removing the casing, disconnect the equipment from the mains.

The manipulation of the internal electronic circuits of the stirrer by unauthorized personnel can cause irreparable damage. Take it to one of the J.P. SELECTA authorized technical services.

CLEANING:

For the cleaning of the different parts, use the following products:

Cleaning of stainless steel: alcohol.

Cleaning of plastics and covers: alcohol with cotton or a soft duster.

SPARE PARTS

To guarantee the safety of the equipment, all spare parts must be purchased from J.P. SELECTA, s.a.u.

DESCRIPTION	Code
Model	7000272
Mantle support	0005086
Mantle support guides	0005087
Control circuit	0038514
Magnet	0015449
Motor	0024283

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by unauthorized personnel by J.P. SELECTA, s.a.u. cancels the guarantee automatically.