FRIGEDOR

3000778

UNIDADES REFRIGERADORAS PARA BAÑOS
IMMERSION COOLERS FOR BATHS

INFORMACIÓN GENERAL**GENERAL INFORMATION**

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The standard equipment consists of the following components:

Pieza	Piece	Código/Code	Cant. / Qty.
FRIGEDOR	FRIGEDOR	3000778	1
Manual de instrucciones	Instruction manual	80016	1

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*** Tanks:**

Thermally insulated to withstand low temperatures. Insulated with injected foam of expanded-polyurethane to avoid condensations between -40°C and +90°C.

Código / Code	Capacidad litros / Capacity litres	Medidas útiles cm. / Tank dimensions cm.			Medidas exteriores cm. / Overall dimensions cm.			Peso Kg. / Weight Kg.
		Alto / Depth	Ancho / Length	Fondo / Width	Alto / Height	Ancho / Length	Fondo / Width	
6003901	9	15	29	22	24	40	34	6
6003921	20	20	48	29	24	61	41	9

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Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Código / Code		3000778
Temperatura °C / Temperature °C		-20 / +20
Medidas exteriores cm. / Overall dimensions cm.	Alto / Height	41
	Ancho / Length	21
	Fondo / Width	34
Potencia frigorífica / Cooling power		a -20°C = 50W
Potencia HP / Power HP		1/5
Consumo W / Consump. W.		285
Peso Kg. / Weight Kg.		14

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EQUIPMENT DESCRIPTION

FRIGEDOR units allow temperatures up to -20°C.

Table-top design with a sealed compressor, a ventilated condenser and a spiral cooling coil in AISI 304 stainless steel.

The quantity of coolant gas is 190gr of R-134a.

COIL DIMENSIONS:

- * Line length: 900 mm.
- * Coil length: 150 mm.
- * Ø coil: 45 mm.

INSTALLATION

Place the machine near a suitable mains connection.

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Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the machine.

Do not use the unit to cool highly inflammable liquids.

Do not use the machine if it is NOT earthed.

If you have to change the plug, bear in mind the following:

Blue cable:	Neutral.
Brown cable:	Phase.
Yellow/green cable:	Earth.

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OPERATION

Turn on the switch with indicator lamp (green).

Place the cooling coil into the tank with liquid which has to be cooled.

If the machine switches off, you must wait for 3 minutes to start it again.

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MAINTENANCE

Before removing the cooler casing, disconnect the machine from the mains.

The guarantee will not have any effect if the machine is manipulated by somebody who has not been expressly authorized by J.P. SELECTA.

CLEANING:

Remove the metallic casing and clean the condenser wings with the help of a brusher or compressed air. It is recommended to do it periodically.

SPARE PARTS

To guarantee the safety of the equipment, all spare parts must be purchased from J.P. SELECTA, s.a.

Código / Code		3000778
Alargo Schuko	<i>Schuko cable</i>	7002
Tirador unidad de frío	<i>Cool unit knob</i>	15490
Interruptor	<i>Switch</i>	20017
Grupo unidad refrigeradora	<i>Refrigerator unit group</i>	24040
Pinzas Jebe	<i>Jebe clamps</i>	27515
Retenedor cable	<i>Cable stopper</i>	28004

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GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, S.A.

Any manipulation of the machine by unauthorized personnel by J.P. SELECTA, S.A. cancels the guarantee automatically.

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