**CENCOM 2**

**7002240**

**CENTRÍFUGA ANGULAR PARA 6 TUBOS DE 15ML**

**ANGULAR CENTRIFUGE FOR 6 X 15 ML TUBES**

## INDICE

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# ENGLISH VERSION

## 1. GENERAL INFORMATION

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.
- 2) Do not install or use the equipment without reading this handbook before.
- 3) This handbook must always be attached to the equipment and it must be available for all users.
- 4) If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta, s.a.u.
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.
- 7) Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.

## 2. PACKING LIST

The standard equipment consist of the following components:

<i>Piece</i>	<i>Code</i>	<i>Qty</i>
CENCOM 2 Centrifuge	7002240	1
Bucket for glass tube		6
Bottom rubber for glass tube		6
Bucket for FALCON tube		6
Bottom rubber for FALCON tube		6
Glass tube 15ml 16.5x100	47105	6
Instruction manual	80044	1

## 3. TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

<i>Max. volume</i>	<i>Max. tube capacity</i>	<i>Overall dimensions</i>	<i>Maximum speed</i>	<i>Maximum R.C.F.</i>	<i>Regulation</i>	<i>Consump. W.</i>	<i>Weight Kg.</i>
90 ml	15 ml	Ø246X240mm	4000 rpm	1790 g	8 positions	92	5

## 4. EQUIPMENT DESCRIPTION

CENCOM 2 is a benchtop centrifuge with an angle head of 6 x15ml tubes. It is supplied with a safety cover and 8 speed positions.

## 5. INSTALLATION

Place the centrifuge on a flat, horizontal, level surface.

Connect the cable into the mains.

**CAUTION!!! IMPORTANT FOR YOUR SAFETY**

**Do not use the equipment to centrifuge liquids which are susceptible to form explosive mixtures.**

**In case of centrifuging microbiological products susceptible to contamination or being contaminated, it is advisable to use tubes with a stopper.**

**Be sure that the voltage supply is the same as the one indicated on the characteristics plate of the equipment.**


**Do not use the equipment if it is NOT EARTHED.**

**If you have to change the plug bear in mind the following:**

<b>Blue cable:</b>	<b>Neutral.</b>
<b>Brown cable:</b>	<b>Phase.</b>
<b>Yellow/green cable:</b>	<b>Earth.</b>

## 6. OPERATION

### CONTROL PANEL:

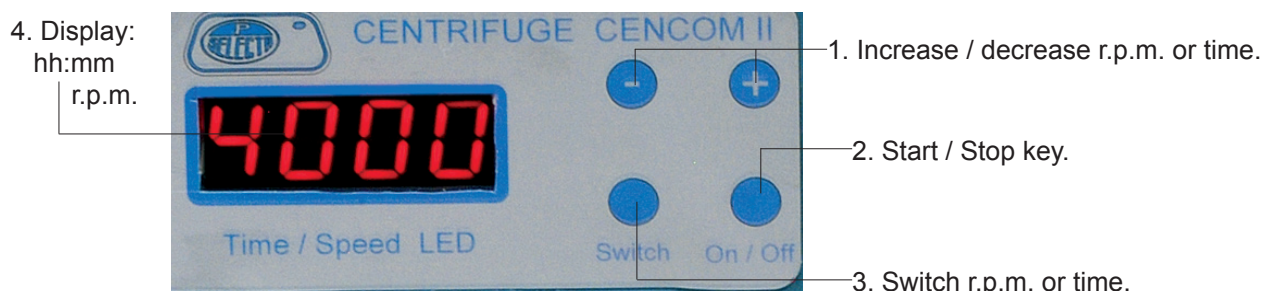## **STARTING UP:**

1. *Connect into mains.*
2. *Put the tubes into the head in a balance way. If all buckets are not used, take notice that the load must be symmetrically balanced.*
3. *Close the cover.*
4. *Select the speed. (r.p.m.):*  
1000, 1500, 2000, 2500, 3000, 3500, 4000.
5. *Select the centrifugation time (hh.mm).*
6. *Press **On / Off** to start centrifugation.*

## **STOP:**

*Press **On / Off** and wait for a few minutes before opening the door. Do not place tubes and remove if the head is not completely stopped.*

*Leave the head to stop by itself. Do not use any tool or your hand to brake the head.*

## **SAFETY:**

*There is a microswitch which disconnects the motor when opening the cover (in case of stop omission of the motor)*

**CAUTION!! DO NOT OPEN THE COVER WHEN THE MOTOR IS OPERATING**

## **7. MAINTENANCE**

***Before removing the cover, disconnect the equipment from the mains.***

### **CLEANING:**


*For the cleaning of the different parts, use the following products:  
Cleaning of casing, plastic and painted metals: Alcohol with cotton duster.*

## **8. GUARANTEE**

*This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.*

*Any manipulation of the equipment by unauthorized personnel by J.P. SELECTA, s.a.u. cancels the guarantee automatically.*

## 9. PLACING TUBES

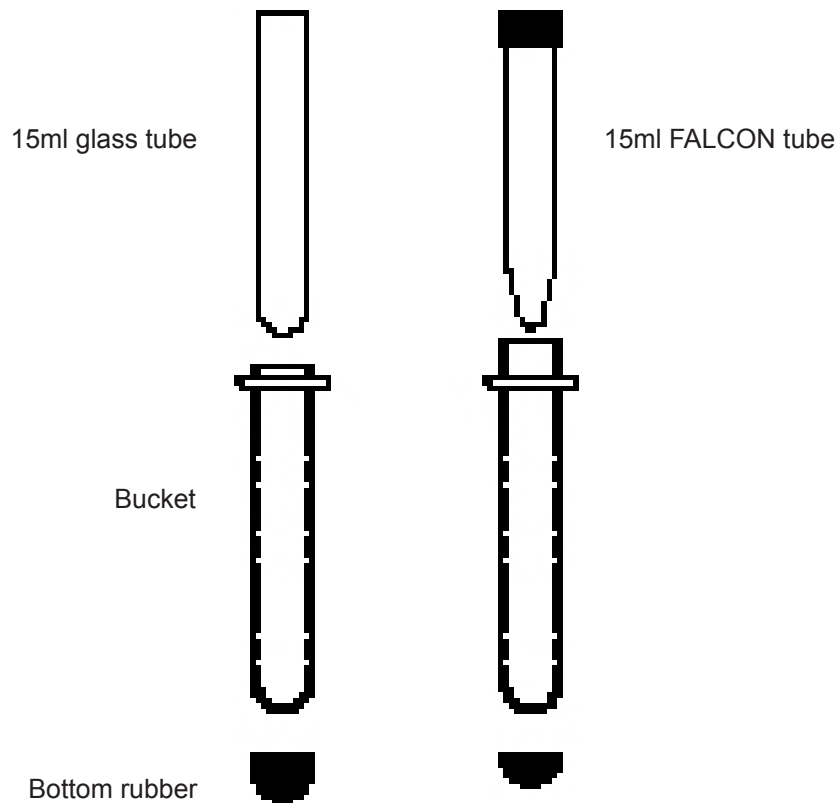## 10. ERRORS TABLE

Malfunction Solutions - Cencom II			
Error code	Code Meaning	Reason	Resolution
E1	Unbalance protection	Unbalanced protection input switch closed	Check imbalance detection switch Unplug the unbalanced protection wire, no alarming occurred
E2	Over speed	Speed measurement interference or speed rise too fast	Check Speed measuring wire or reduce the speed up
E3	Door cover protection	Door cover input signal wire is not connected properly or Door cover switch fault	Check the door cover input signal wire and switch
E6	Over current	Current over rated value	Speed up too fast under heavy load, reduce speed increasing Drive board failure External voltage too high
E7	No speed measurement	No speed signal input The speed sensor damaged or wire connection loose	Check speed sensor and make sure the wire connection is good There are many reasons for the input of no speed measurement signal Analyze the cause separately
E11	The set speed can't be reached	Insufficient motor power External voltage too low	Ensure load and control system power match Ensure the power supply is normal
	No speed measurement of motor after starting, please turn the motor manually. Observe whether the LED beside the speed measuring interface flashes, If it does not flash, the speed sensor may be damaged. If flashing, contact after-sales service.		