**AGIMATIC - HS**  
**AGIMATIC - HL**

**7001005**  
**7001006**

**AGITADORES MAGNÉTICOS POTENTES**  
**POWERFUL MAGNETIC STIRRERS**

## ÍNDICE

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# ENGLISH VERSION

## GENERAL INFORMATION

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.
- 2) Do not install or use the equipment without reading this handbook before.
- 3) This handbook must always be attached to the equipment and it must be available for all users.
- 4) If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 2009/104/CE , the manufacturer is not responsible for the damage that can occur.
- 7) Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.

## PACKING LIST

The standard equipment consist of the following components:

Piece	Code	Quantity
AGIMATIC - HS	7001005	1
Magnet	1001063	1
AGIMATIC - HL	7001006	1
Magnet	1001064	1
Instructions manual	80109	1

## ACCESSORIES

### MAGNETIC STIRRING BARS

Code	Dimensions
1000009	6Ø x 10mm
1000012	4Ø x 12mm
1000019	7Ø x 20mm
1000032	8Ø x 32mm
1000039	8Ø x 40mm
1000045	10Ø x 45mm
1001061*	12Ø x 75mm
1001062*	16Ø x 102mm
1001063*	16Ø x 127mm
1001064*	19Ø x 150mm

\*For big volumes, with a movable pivot ring.

## TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

MAGNETIC STIRRER	Code	Speed regulation	Stirring power	Overall dimensions cm			Consump.	Weight
				Height	Length	Width		
AGIMATIC - HS	7001005	50 - 980 r.p.m.	Up to 25 litres	12	28	34	100 W	8 Kg
AGIMATIC - HL	7001006		Up to 60 litres	14	44	38	190 W	12 Kg

## EQUIPMENT DESCRIPTION

Magnetic stirrers for volumes from 25 up to 60 litres of low viscosity, depending on model.

With electronic speed regulation and constant torque which is maintained regardless of fluid's viscosity.

## INSTALLATION

Place the apparatus on a flat horizontal, level surface, leaving a space at the back and sides of the apparatus of about 10cm.

### **CAUTION!!! IMPORTANT FOR YOUR SAFETY**

Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the apparatus.


Do not use the apparatus if it is NOT earthed.

If you have to change the plug bear in mind the following:

**Blue cable: Neutral.**  
**Brown cable: Phase.**  
**Yellow/green cable: Earth.**

## OPERATION

### CONTROL PANEL:

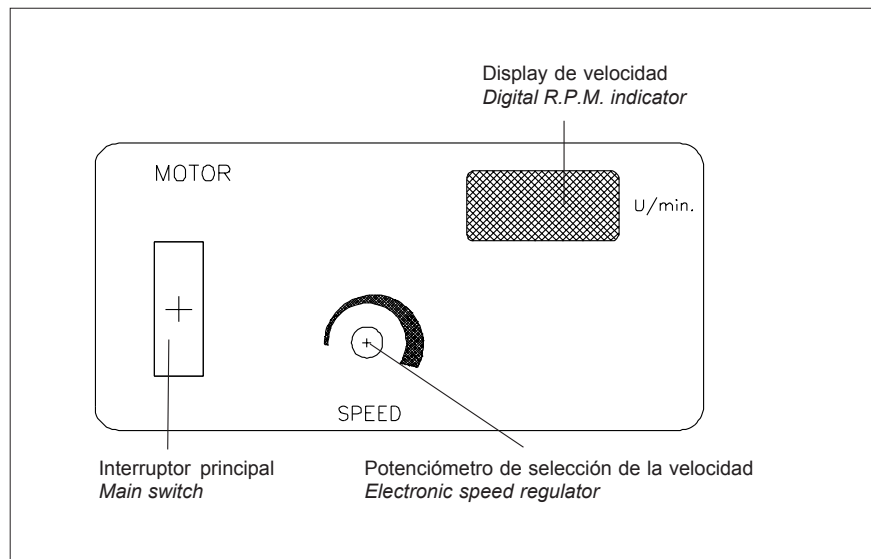- Place the recipe with the adequate magnet in the middle of the top plate.
- Turn the speed regulator at minimum power (anticlockwise).
- Press the main switch in the ON position.
- Turn slowly the speed regulator clockwise until the desired speed.

## MAINTENANCE

Before removing the cover, disconnect the equipment from the mains.

The manipulation of the internal electronic circuits of the equipment by unauthorized personnel can cause irreparable damage. Take it to one of the J.P.SELECTA's authorized technical services.

## SPARE PARTS

To guarantee the safety of the equipment, the spare parts must be purchased from J.P.SELECTA, s.a.u.

Description	Code
Control P.C.B.	38528
Tachometer P.C.B.	29013
Motor	24068

## GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damages caused by incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by unauthorized personnel by J.P. SELECTA, s.a.u. cancels the guarantee automatically.

### Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.