PHARMALOW - S
PHARMALOW - L

2101274
2101275

FRIGORÍFICO INDUSTRIAL
COMMERCIAL REFRIGERATOR

ADVERTENCIAS GENERALES

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SAFETY INSTRUCTIONS

- 1) *To prevent injury or damage to the unit, the appliance should be unpacked and set up by two people.*
- 2) *In the event that the appliance is damaged on delivery, contact the supplier immediately before connecting to the mains.*
- 3) *To guarantee safe operation, ensure that the appliance is set up and connected as described in these operating instructions.*
- 4) *Disconnect the appliance from the mains if any fault occurs. Pull out the plug.*
- 5) *When disconnecting the appliance, pull on the plug, not on the cable.*
- 6) ***Any repairs and work on the appliance should only be carried out by the customer service department, as unauthorised work could prove highly dangerous for the user. The same applies to changing the mains power cable.***
- 7) *Do not stand on the plinth, drawers or doors or use them to support anything else.*
- 8) ***If you have a lockable appliance, do not keep the key near the appliance or within reach of children.***
- 9) *Do not store explosives or sprays using combustible propellants such as butane, propane, pentane etc. in the appliance. Electrical components might cause leaking gas to ignite. You may identify such sprays by the printed contents or a flame symbol.*

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SETTING UP

- 1) *Fit the handle into the attachment holes. By fitting the handle yourself you will help to save on packaging material.*
- 2) ***Avoid positioning the appliance in direct sunlight or near cookers, radiators and similar sources of heat.***
- 3) *The floor on which the appliance stands should be horizontal and level. Compensate for uneven floors with the adjustable feet (SW 22). The appliance can be placed with its back flush to the wall.*
- 4) *Always ensure that there is good ventilation and that the outward flowing air is able to escape.*

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CONNECTING TO THE MAINS

Power supply (a.c.) and voltage at the operating point must comply with the details on the type plate, which is located on the outside of the appliance. The socket must be fused with a 10 A fuse or higher.

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To ensure trouble-free operation, do not connect the appliance to the power supply until about 6 to 8 hours after delivery. Connect the appliance through a properly earthed plug and socket.

The socket should be easily accessible and not directly behind the appliance.

The wires in the mains lead are coloured in accordance with the following code: green/yellow = earth, blue = neutral, brown = live.

Warning! This appliance must be earthed.

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SWITCHING THE APPLIANCE ON AND OFF

You are recommended to clean the appliance before switching it on for the first time (see "Cleaning"). When it is connected up (plugged into an approved wall socket) the appliance is ready for operation.

To switch the appliance off: pull out the mains plug

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SETTING THE TEMPERATURE

Using a coin, turn the arrow on the temperature control to a setting between "1" and "7".

Setting "1" = lowest cooling setting, highest temperature.

Setting "7" = highest cooling setting, lowest temperature.

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COOLING

This appliance can be used for chilling all types of materials. The grid shelves can be moved to accommodate different height packages or objects.

To remove the shelves the door must be opened wider than 90°.

Substances which give off or absorb odours and flavours as well as liquids should always be stored covered or in closed containers. High-percentage alcohols should be tightly sealed and stored upright.

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The ventilation slots on appliances with recirculated air fans on the inside must be not covered.

NB: This commercial appliance is suitable for almost universal use.

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SHUTTING YOUR APPLIANCE DOWN

If your appliance is to be shut down for any length of time, switch it off and disconnect the plug. Clean the appliance and leave the door open in order to prevent un pleasant smells.

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DEFROSTING

The refrigerator compartment defrosts automatically. The water that forms on the rearwall drains into a reservoir at the back of the appliance and evaporates automatically through the compressor heat. Ensure that the water can flow without hindrance through the drain hole (see "Cleaning").

A layer of frost or ice can form on the cooling elements when the appliance has been operating at the coldest setting for a long time.

*If a **layer of ice** has formed, the cooling elements should be defrosted and the inside surfaces cleaned.*

Do not use any mechanical devices or other artificial aids for defrosting other than those recommended by the manufacturer.

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CLEANING

Before cleaning always switch off the appliance. Pull out the mains plug.

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1. Clean the inside, equipment parts and outer walls with lukewarm water and a little detergent.

Do not use steam cleaners because of the risk of injury and damage.

Do not use abrasive or acid cleaners or chemical solvents. We recommend the use of an all-purpose cleaner with neutral pH.

2. Ensure that no cleaning water penetrates into the electrical components or ventilate grille.

3. Dry all parts well with a cloth.

4. The dust should be removed from the refrigeration unit and heat exchanger-metal grid at the back of the appliance - once a year.

5. Clean the drain hole in the refrigerator compartment with a thin object, e.g. cotton swab or the like.

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MALFUNCTIONS

Your appliance is designed and manufactured for a lengthy, reliable operating life with no malfunctions. If a malfunction nonetheless occurs during operation, check whether it is due to an operating error. Please note that even during the warranty period the resultant servicing costs in this case will have to be borne by the owner.

You may be able to rectify the following faults by checking the possible causes yourself:

1. Appliance does not function:

- *Is the appliance switched on?*

- *Is the plug correctly fitted in the mains socket?*

- *Is the socketed fuse intact?*

2. Loud running noise:

- *Is the appliance set up firmly on the floor?*

- *Does the appliance cause nearby items of furniture or objects to vibrate?*

Ensure that the appliance is stable and that all parts at the back can vibrate freely.

Please note that noises caused by the refrigerant circuit cannot be avoided.

3. The temperature is not low enough:

- *Is the temperature setting correct (see "Setting the temperature")?*

- *Does the separately installed thermometer show the correct reading?*

- *Is the ventilation system working properly?*

- *Is the appliance set up too close to a heat source?*

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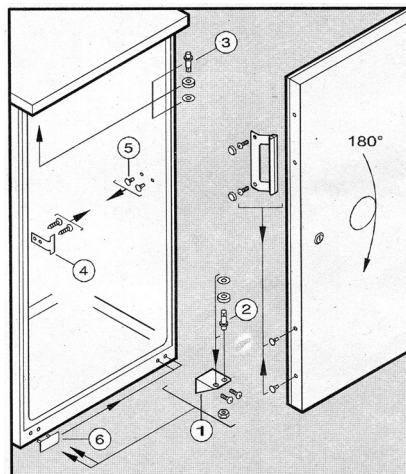
3. *The temperature is not low enough:*
- *Is the temperature setting correct (see "Setting the temperature")?*
 - *Does the separately installed thermometer show the correct reading?*
 - *Is the ventilation system working properly?*
 - *Is the appliance set up too close to a heat source?*

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1. *Close the door and unscrew the hinge bracket ① (support the door if necessary).*
2. *Retire the door downwards.*
3. *Unscrew the hinge pin ② from the hinge bracket ① and screw into the adjacent retaining hole.*
4. *Unscrew hinge pin ③ and screw in on opposite side.*
5. *Attach the closure ④ and plugs ⑤ on the opposite side.*
6. *Rotate door by 180°.*
7. *Screw the handle onto the door using the holes provided.*
8. *Lift off the cover ⑥ and seal the exposed retaining holes on the other side with it.*
9. *Mount door on hinge pin ③ and close.*
10. *Insert hinge bracket ① in lower door mounting and screw onto plinth.*
11. *If the door is not flush with the body of the appliance, adjust it using the slots on the hinge bracket ① (ensuring that all screws are securely fastened).*

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