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The standard equipment consists of the following components:

<b>Descripción</b>	<b>Cant.</b>	<b>Description</b>	<b>Qty.</b>
- Vidrio condensador	1	- <i>Condenser</i>	1
- Tapa de vidrio	1	- <i>Glass lid</i>	1
- Calderín inox	1	- <i>Vessel</i>	1
- Junta de la tapa	1	- <i>Lid's ring</i>	1
- Pinza cono 34/35	1	- <i>Stopper / "O" ring gasket</i>	1
- Manguera de silicona entrada de agua	1	- <i>Feed tube</i>	1
- Manguera de silicona salida desagüe	1	- <i>Drain tube</i>	1
- Manguera de silicona salida agua destilada	1	- <i>Distillate tube</i>	1
- Manguera de silicona nivel	1	- <i>Level tube</i>	1

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## GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please notify to your J.P. Selecta supplier immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubt or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lack of maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the equipment liquids which can give off vapours capable of making explosive mixtures.*

## GUARANTEE

*This product is guaranteed for one year. Guarantee does not cover damages caused by an improper use or by causes beyond J.P. SELECTA, S.A. control.*

*Any manipulation of equipment made by personnel not authorized by J.P. SELECTA, S.A. automatically cancels the guarantee.*

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**ATTENTION!: This equipment do not have fuse. A circuit breaker and a differential switch must be installed at mains.**

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**ATTENTION**

Since distilled water is in contact with the atmosphere before the conductivity measurement can be carried out, its value is actually much less (up to 10 times less) than what it is actually measured. The measurement of any ultrapure water in contact with atmosphere has a conductivity of 1-2 µS/cm and a pH of 5, due to a small amount of CO<sub>2</sub> (0.05 ppm) absorbed from the atmosphere.

	Conductivity (uS/cm)	Resistivity (MOhm · cm)	Silicates (mg/l)	Heavy metals (mg/l)	Reduced Permanganate (min)	Sodium (mg/l)	Calcium hardness (mg/l)	Ammonium (mg/l)	Bacteria (u.f.c./ml)	pH (at 25 °C)
Once distilled	10 – 2	0,1 – 0,5	1 – 0,5	1 – 0,5	30	5 – 2	3 – 1	0,01	<10	5 – 6,5
Bidistilled	2 – 1	0,5 – 1	0,7 – 0,1	0,8 – 0,1	30	1 – 0,5	0,3 – 0,1	0,01	<10	5 – 6,5

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*To install Aquasel distilling you must follow the next items (see picture):*

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1. *Put the lid's ring supplied in the groove of the vessel.*

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2. *Place the glass lid on the ring and lock it with the clips.*

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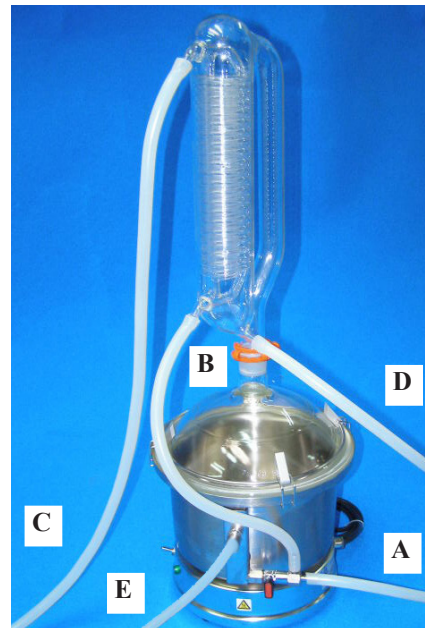
4. *Connect pipes as showed in the pictures.*

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
- A. Distillation and refrigeration water feeding.  
 B. Refrigeration water inlet  
 C. Refrigeration water outlet  
 D. Distilled water outlet  
 E. Overfeeding water outlet*

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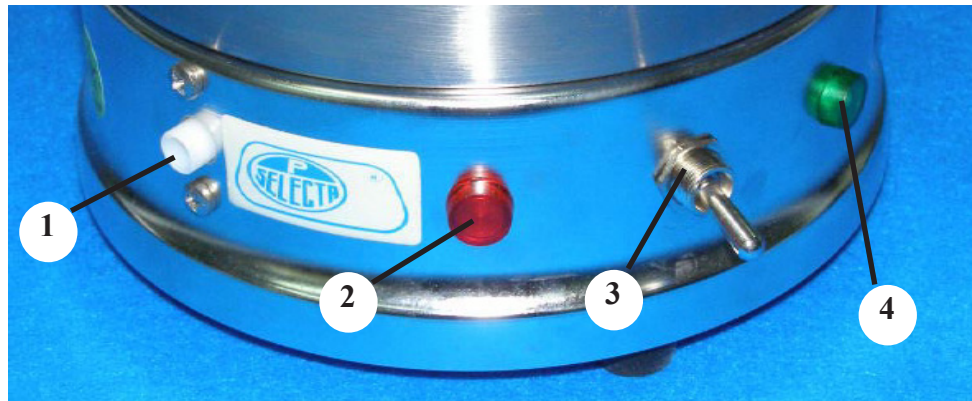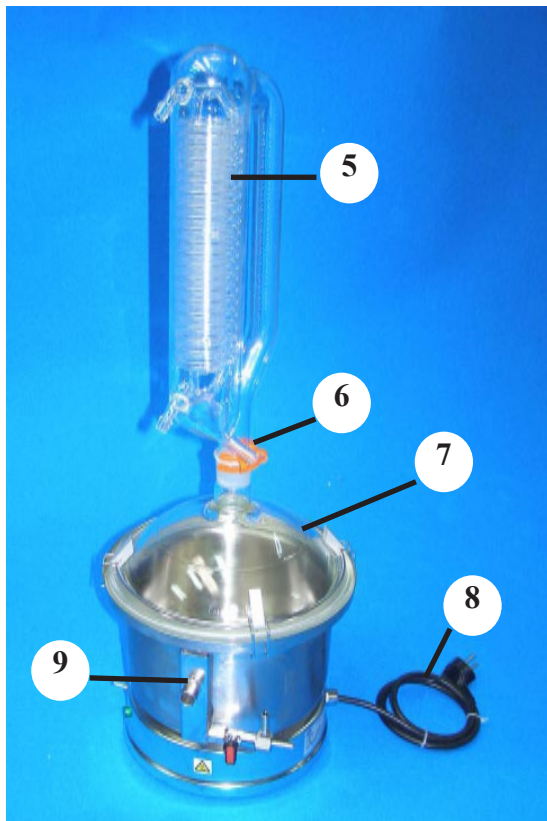
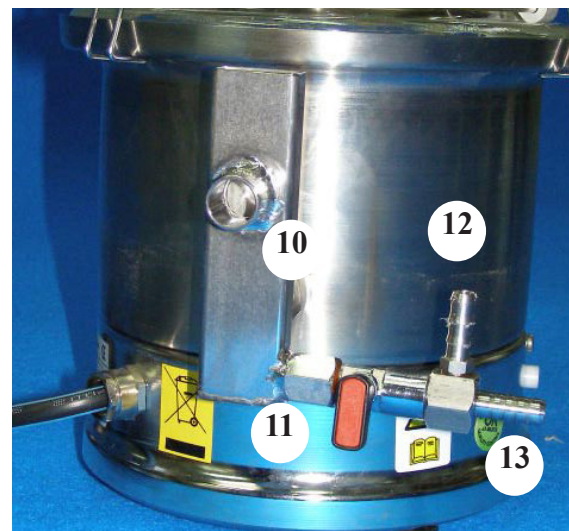
Detalle nivel constante  
Constant level detail

5. Glass condensor
6. Looking condenser grip
7. Bath glass lid
8. Power supply cord
9. Constant level
10. Overfeeding water outlet
11. Water level regulation valve
12. Cooling water
13. Water inlet

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## STARTING UP

1. *Keep in mind that the installation was done correctly.*
2. *Turn on the water tap until the water covers the heater elements of the bath.*
3. *Turn off the tap.*
4. *Turn on the main switch (3). The green pilot lamp (4) will light up.*
5. *Keep the level regulation valve (11) turned off.*
6. *When the water begins to boil, turn on the tap.*
7. *Turn on slowly the regulations valve (11) until the water outfalls by the hose (E) with a very small outflow.*  
*Keep in mind that the flow regulated by the valve (11) only must be enough to back fill the water evaporated by the distillation.*

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**IMPORTANT: Be sure that the heater elements are submerged before connect it to the mains.**

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## MAINTENANCE

*The calcium depositions on the heater elements lose calorific power.*  
*Depending of the water's hardness, the lime must be cleaned frequently. To remove the lime, one normally uses acetic acid to 20% or a ready-made product.*

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**IMPORTANT: Before any operation disconnect the heater elements from de mains.**

## RECAMBIOS

## SPARE PARTS

Descripción Description	Códigos Codes	
	4903004	4903006
Resistencia calefactora Heater element	39232	39229/39229
Junta calefactora Lid's ring	21155	21156
Clip unión 34/35 y pinza Hoffman O-ring gasket and stopper		1001100/1000064
Tapa vidrio Glass lid	47025	47027
Condensador de vidrio Glass condenser	47026	47028