ULTRASONS-UB-1488

3001488

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***ULTRASONIC CLEANERS
WITH ELEVATOR AND AUTOMATIC DRYER***

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please notify to your J.P. Selecta supplier immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubt or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lack of maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

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PACKING LIST

The standard equipment consist of the following components:

- Ultrasonic cleaner Cod. 3001488
- Basket Cod. 09449
- Wire Cod. 07001
- Instruction manual Cod. 80166

EQUIPMENT DESCRIPTION

Ultrasonic cleaners with basket elevator and automatic drying at the end of cycle.

The principle of ultrasonic cleaning consist on the use of high frequency sound waves (40 KHz) produced by a generator through a transducer, which propagates them mechanically inside the tank, producing a cavitation effect which leads to the formation of millions of low pressure microscopic bubbles which carry out a molecular cleaning.

Parts containing sites with complex acces can be cleaned without disassembly because the cavitation effect reach wherever the cleaning solution penetrates.

TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

| | | |
|--|--------------|--------|
| Capacitat (litros)/Capacity (litres) | | 6 |
| Medidas útiles (cm) Tank dimensions (cm) | Alto/Height | 110 |
| | Ancho/Length | 265 |
| | Fondo/Width | 125 |
| Medidas exteriores (cm) Overall dimensions (cm) | Alto/Height | 325 |
| | Ancho/Length | 440 |
| | Fondo/Width | 340 |
| Peso (kg)/Weight (kg) | | 15 |
| Potencia generador (W)/Generator power (W) | | 150 |
| Potencia secador (W)/Dryer power (W) | | 1000 |
| Limpieza temporizada (min)/Clean timer (min) | | 15 |
| Secado temporizado (min)/Drying timer (min) | | 15 |
| Filtro de red/Mains filter | | Sí/Yes |

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INSTALLATION

Place the apparatus on a flat, horizontal, level surface, trying to leave a space at the back of the apparatus of 10 cm.

To make easier package and transport the apparatus is delivered with elevator in lower position (cleaning position). When cleaner is connected at mains and the main switch is switched on, the elevator up automatically.

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Do not use the cleaner with inflammable liquids which can give off vapours capable of making explosive mixtures.

Be sure that the voltage supply is the same as the one indicated on the characteristics plate.

Do not use the apparatus if it is not earthed.

If you change the plug bear the following in mind:

Blue cable: Neutral.
Brown cable: Phase.
Yellow/green cable: Earth.

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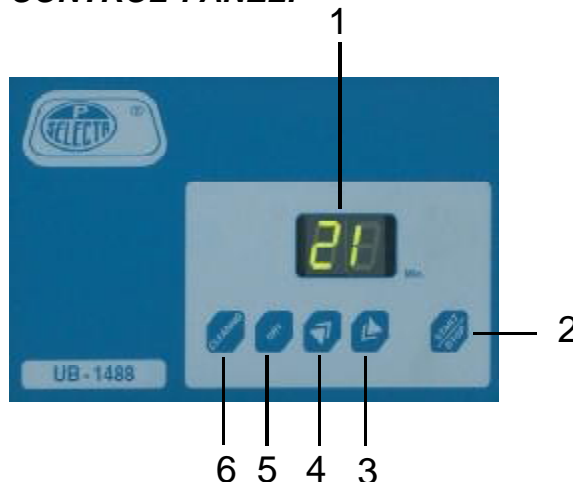
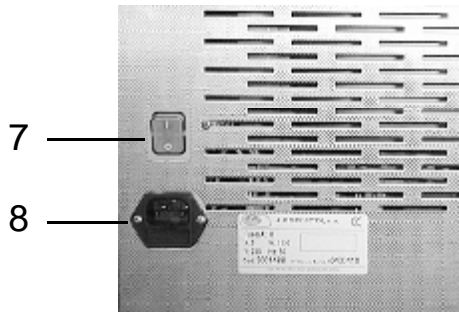
1. *Switch on the mains switch (7) on the rear part*
2. *Set cleaning time pusching CLEANING (6) and at the same time, (3) and (4) keys to increase or decrease time in minutes. Highest time 15 minutes.*
3. *Set drying time pusching DRY (5) and at the same time, (3) and (4) keys to increase or decrease time in minutes. Set 0 minutes to avoit drying cycle. Highest time 15 minutes.*
4. *Put the material into the basked being in mind that weight of load be lower than 2,5 kg.*
5. *Push START/STOP key (2) to start clean process and dry.*

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Clean cycle begins submerging the basket in cleaner solution.
When clean process ends basket emerge and if drying time is set drying cycli starts.
When all cycles are finished it sounds and acoustic signal

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- * *Do not overload the cleaner. Avoid placing heavy instruments directly on the bottom of tank.*
- * *Clean process produces heat so clean solution gets warm.*
- * *The first it has to clean fragile or valuable pieces, you should try to clean only one of them before continue cleaning the others. Never use ultrasonic cleaners with organic pearls.*

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CLEANING SOLUTIONS:

Among cleaning solutions we can find two main kinds: those which have water as the basic element in which a detergent is added, and other solvents. The detergents which are mixed with water may be solid or liquid. Water is the most universal cleaning type, as it is able to dissolve many organic substances. Although water may not dissolve such organic substances as oil, grease or wax, but it can be used to carry the cleaning detergent that act on these products.

Water also acts as environment of dispersion for organic substances which cannot be dissolved but which may be carried by its.

Most of cases, and mainly when there are no large traces of grease or oils, a standard domestic liquid detergent may be proper.

When volume and type of dirt is such that aqueous solutions do not attack it at all or do it very slowly (wax, oils or greases), specific solvents have to be used.

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* **Detergent BIO-SEL:** Specially for cleaning with **ULTRASONS** ultrasonic cleaners. We advise of this special detergent by its excellent cleaning power, it is biodegradable and free of phosphates. It is very useful for cleaning electronic components, pieces of glass, metal, plastic, rubber, jewels and materials with hard surfaces in general.

For cleaning long objects whose length may exceed that of the tank, such as tubes, pipettes, etc., it has been proved that they are submitted to the same cavitation effect by introducing them in a test tube filled with liquid and placed vertically in the cleaner.

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
EMPTYING TANK:

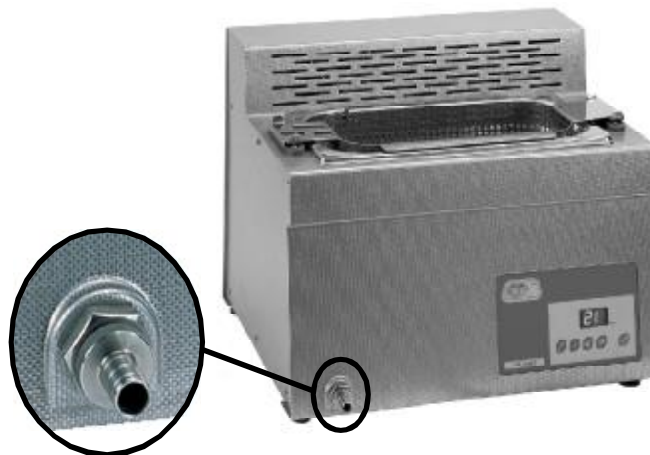To empty the tank it is recommended to connect a hose at the outlet placed on the front part of cleaner and unscrew the nut to turn the emptying valve on. The tank content can be collected in a recipient in order to treat the product properly or be emptied in waste pipe.

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MAINTENANCE

Before removing the side covers disconnect the apparatus from the mains.

The guarantee will not have any effect if the apparatus is manipulated by somebody who has not been expressly authorized by J.P.SELECTA.

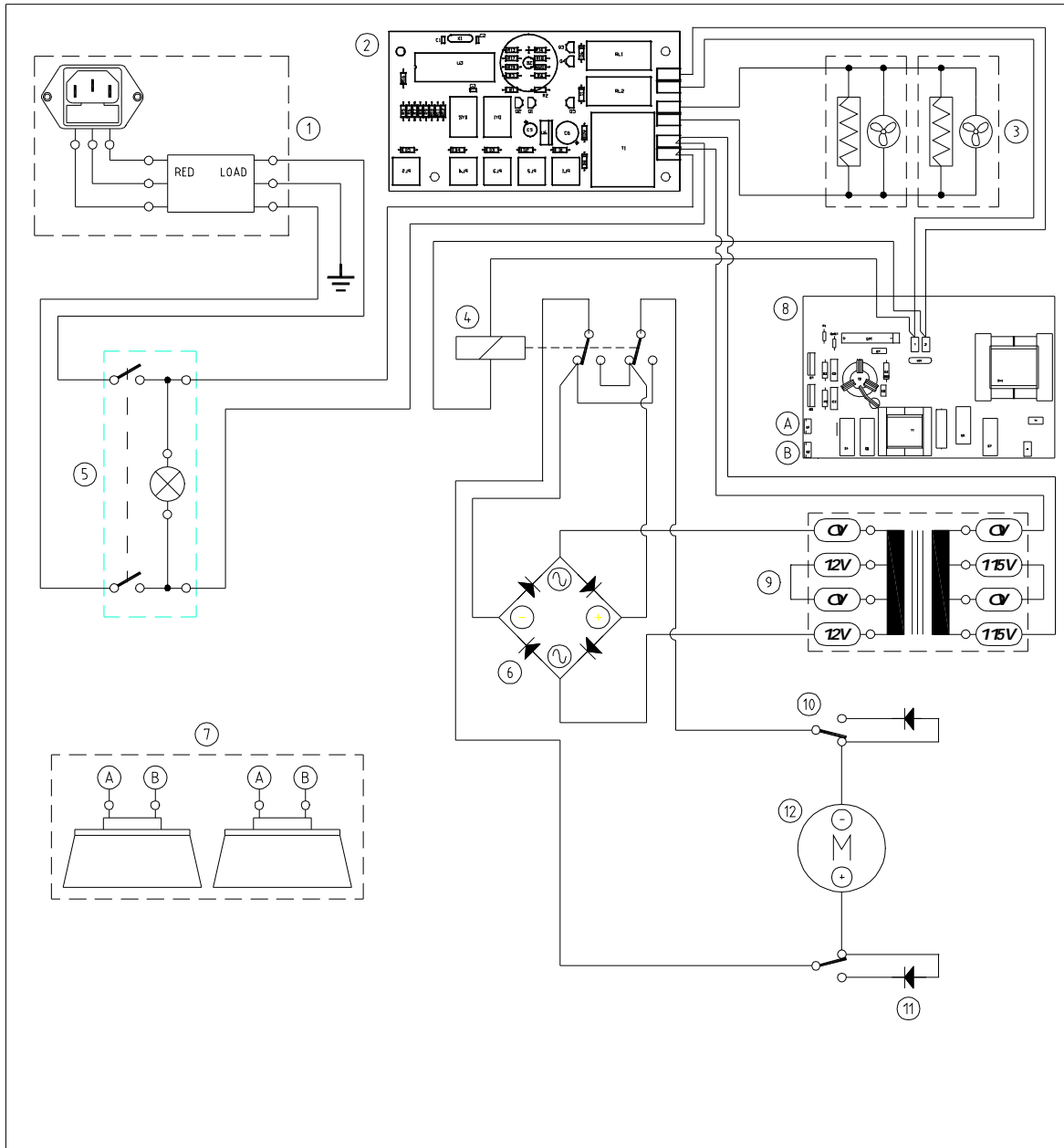
CLEANING:

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol.

Cleaning of plastics and covers: Alcohol with cotton or a soft duster.

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GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use of causes beyond the control of J.P. SELECTA, S.A..

Any manipulation of the apparatus by unauthorized personnel by J.P. SELECTA, S.A. cancels the guarantee automatically.

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