4100050 Sealcom-V

Selladora de bolsas al vacío

Sealing bags under vacuum

Indice

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
GENERAL INFORMATION

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.
- 2) Do not install or use the equipment without reading this handbook before.
- 3) This handbook must always be attached to the equipment and it must be available for all users.
- 4) If you have any doubts or enquiries, please contact with your supplier or with J.P. Selecta's technical service.
- 5) **IMPORTANT! J.P. Selecta, s.a.u. WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.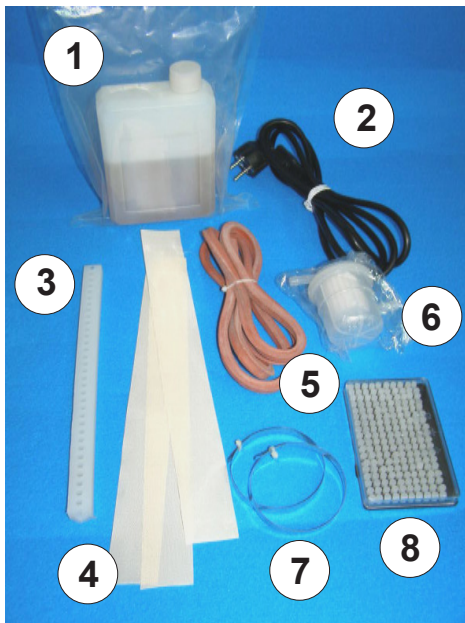

PACKING LIST

The standard equipment consists of the following components:

Sealing machine	4100050
Vacuum pump oil (spare part)	1
Power supply wire	2
Silicone shoe (spare part)	3
Ignifuge heater cover (spare part)	4
Lid gasket (spare part)	5
Air pump filter (spare part)	6
Heating elements (spare part)	7
Letterpress set	8
Instruction manual	80212

TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Chamber dimensions:

Length	32 cm
Width	28 cm
Height	05 cm

Overall dimensions:

Height	28 cm
Width	32 cm
Depth	52 cm

Weight:	35 kg
Heating power:	370 W
Voltage:	230 V
Intensity:	0.39 A
Vacuum level:	-0.1 MPa
Sealing width:	10 mm
Sealing max. length:	260mm

EQUIPMENT DESCRIPTION

The bag seling Vacuum Packing is a device easy to use and adaptable to different uses because it allows automate vaccum time, sealing time, cooling time, 3 hetting levels and cool sealing.

Bag stamp ability

INSTALLATION

Place the sealing machine on a horizontal and level surface keeping in mind the weight of the device: 36 kg

Before to start up the device check the oil level of the pump as well is indicated on the section "Starting up".

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Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the apparatus.

Do not use the equipment if it is NOT earthed.

If you have to change the plug keep in mind the following:

Blue cable: Neutral.




Brown cable: Phase.

Yellow/green cable: Earth.

OPERATION

CONTROL PANEL AND OTHER ELEMENTS:

- 9 Main switch
- 10 Power supply socket
- 11 Oil level vacuum pump display
- 12 Pump oil inlet
- 13 Air suction holes
- 14 Emergency stop key 
- 15 Time set key 
- 16 Increase time set 
- 17 Decrease time key 
- 18 Heating level set key 
- 19 Vacuum run led
- 20 Sealing run led
- 21 Cooling indicator led
- 22 Working indicator led
- 23 Display
- 24 Heating level indicators
- 25 Heating element

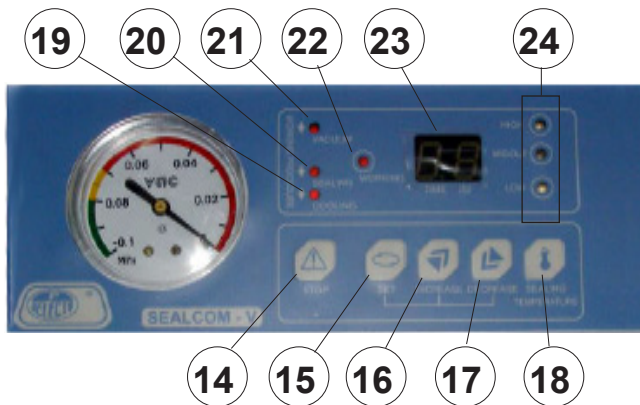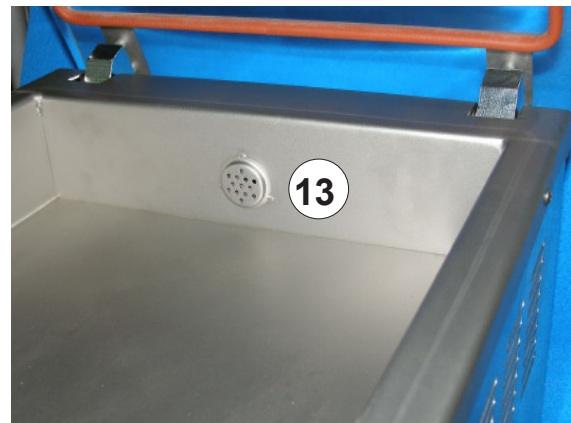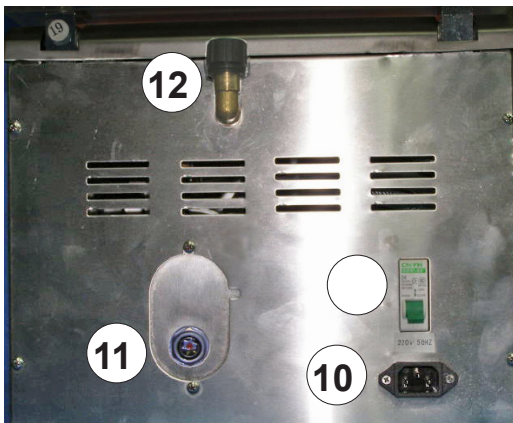STARTING UP:

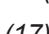
THE PUMP CANNOT WORK IF THE PUMP DOES NOT HAVE A SUITABLE OIL LEVEL

- * The maximum level is the 3/4 of the round window
- * The minimum level is the bottom of the red point

The working parameters must be set before starting the sealing process with the methacrylate lid opened.

- 1 Turn on the main switch (9) on the rear part. The display (23) shows “__”. The sealing machine is “stand by”. The lid must remain opened, if not, it will start the vacuum and sealing process.
- 2 Press the  key (15) to set the vacuum time (between 0 and 99 seconds) using the  (16) and  (17) keys. The VACUUM (21) led will light up.
- 3 Press the  key (15) again to save the value and follow with the next step. The SEALING led (20) will light up.
- 4 Set the sealing time (between 0 and 9.9 seconds) using the keys  (16) and  (17).
- 5 Press the key  (15) again to save the value and follow with the next step. The COOLING led (19) will light up.

- 6 Set the cooling time (between 0 and 9.9 seconds) using the keys  (16) and  (17).
- 7 Press the  (15) once again to exit of the set mode. The display (23) will show "Ed".
- 8 Set the heating level:
Three heating levels and cool sealing can be set. Set it by using the  key (18). The leds HIGH (high level), MIDDLE (middle level) or LOW (low level) will light up. If none of the temperature levels (24) lights up, it means cool sealing.

**TO STOP THE PROCESS
USE THE  (14) AS EMERGENCY STOP**

WORKING:

Once the working parameters are set, the work can be started:

- 1 Place the bag containing the product with the opened end on the heating element and hold it with the transverse bar.
- 2 Close the lid by locking it with the transverse bolt and the cycle will automatically start up with the set parameters: vacuum time, sealing time, cooling time and heating level.
- 3 When the cycle ends, the chamber is automatically pressurized. If the chamber was not locked with the bolt, it opens itself.

MAINTENANCE

Before removing the cover, disconnect the equipment from the mains.

The control elements and the heating elements must only be manipulated by qualified personnel.

CLEANING:

It is advisable to clean the vacuum chamber with a cloth damped with soapy water. Rinse carefully using the same cloth.

Keep the air suction holes cleaned and clear.

Before starting up check, the pump oil level as well is indicated on the section "Starting up". If the level is too low, fill in with the advised oil through the inlet (12)

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. Selecta, s.a.u. Any manipulation of the equipment by unauthorized personnel by J.P. Selecta, s.a.u. cancels the guarantee automatically.