CENTRIFUGAS DE CONTROL ELECTRONICO-  
DIGITAL  
Y MOTOR DE INDUCCION  
ELECTRÓNICO-DIGITAL CONTROL CENTRIFUGES  
WITH INDUCTION MOTOR

CENTROLIT II-BL 7002357

## Índice

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English version

# 1 Safety

The equipment comes complete with safety features.

This manual indicates areas of possible risk.

## 1.1 Safety Icons

The safety icons identify risk situations and safety measures that should be adhered to.

The icons make reference within the paragraph, marked as a grey line.

### Risk danger

Risk danger

Beware, follow the instructions as indicated.

### Electrical risk

Risk of electrical shock if accessing zones are shown with this icon.

Beware, follow the instructions as indicated.

### Important information

- Important information in obtaining the best results or optimum performance of the equipment.
- Important information to extend the life of the equipment and maintain its optimum performance.

## 1.2 Risks that the operator can be subjected to:


- Blows risk.
- Electrical risk.

## 1.3 Qualified users

This equipment should only be used by personnel who are qualified users.

This equipment should only be used by personnel who have understood this manual or have previously adequate instruction of the use of steam sterilisers.

## 2 General information

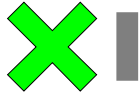- 2.1 Handle the equipment with care. Unpack and check that all items coincide with the delivery note. If you see any discrepancy or damage of any kind, notify your distributor as soon as possible.
- 2.2 Read this manual before using the equipment.
- 2.3 These instructions are an important part of the equipment and should be made available to all users.
- 2.4 If you have any doubts or need to clarity on the use of this instrument, contact your local distributor or the service department of J.P. Selecta, s.a.
- 2.5 **ATTENTION, EQUIPMENT NOT CLEAN AND DISINFECTED SHALL NOT BE REPAIRED.**
- 2.6 Any unauthorised modification, lack of maintenance or deliberate damage to the machine, transgresses the directive 89/655/CEE where the manufacturer will not be responsible for damages that could occur due to undue care and attention.
- 2.7 Do not use the equipment for fluids that can produce vapours which can form inflammable or explosive mixtures.

## 3 Technical specifications.

**Voltage supply 115-230V 50/60Hz according to the characteristics plate indications.**

- Code: 7002357
- Dimensiones (cm):
  - High: 20
  - Width: 28
  - Depth: 34
- Power (W): 170
- Speed max. (rpm): 12000
- RCF max.: 14103 (Depending the head)
- Weight (kg): 6

## 4 Contents list

### 4.1 Standard contents

The standard equipment consists of the following items:

- Centrifuge: 7002357
- Power supply cord 7140
- Instruction manual 80225

### 4.2 Accesorios

Interchangeable heads:

| Head code | For test        | Maximum volume ml. | Tube capacity ml.                     | R.p.m. máx. | R.C.F. máx. (xg) |
|-----------|-----------------|--------------------|---------------------------------------|-------------|------------------|
| 7001704   | Microhematocrit | --                 | 24 of Ø1.4 x 75mm                     | 12000       | 13143            |
| 7001703   | Eppendorf       | 40'5               | 18 of 0.25/0.5ml +<br>18 of 1.5/2.2ml | 12000       | 14103            |

\*MICROHEMATOCRIT READER

Suitable al head 7001704. Direct reading. Code: 7000929

## 5 Introduction

The centrifuge Centrolit II-BL is a tabletop centrifuge with a free maintenance induction motor and electronic digital control of speed and working time, controlled acceleration and deceleration, temporized mode from 1 minute to 60 minutes and safety electromechanic lock.

## 6 Equipment description


### 6.1 Equipment

The Centrolit II-BL centrifuges have the following features:

- Electronic digital control circuit board
- Induction motor (brushless)
- Electromagnetic safety lock

### 6.2 Features


- Regulable speed until 12000 rpm
- Max RCF 14103 g (depending on the head)
- Speed reading on the display (resolution 10 rpm)
- Working time reading on display (resolution 1 min)
- Instant start/stop key to the set speed
- Electronic lid opening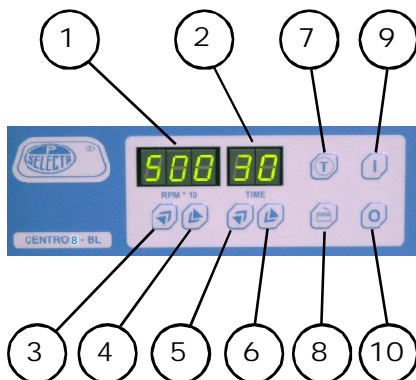

### 6.3 Control panel

1. Speed display (rpm)
2. Time display (min)
3. Increase speed value key
4. Decrease speed value key
5. Increase time value key
6. Decrease time value key
7. Instant start/stop key
8. Start key
9. Stop key

### 6.4 Rear part

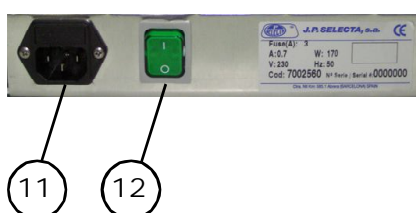10. Power supply socket
11. Main switch

## 7 Installation

### 7.1 PLACEMENT

Place the piece of equipment near of a socket keeping in mind its amperage on a leveled surface enough resistant to support the weight indicated on the table of the section 3 (Technical features), and leaving an space of 10 cm by the rear part and both laterals. If it is a table make sure that its feet are able to absorb the centrifuge vibrations.

DO NOT INSTALL THE CENTRIFUGE IN AREAS WHERE INFLAMMABLE LIQUIDS ARE STORED OR SPECIAL PROTECTION AREAS.

### 7.2 POWER SUPPLY CONEXION

Make sure that the voltage given by the socket is the indicated on the feature label attached on the rear part .

Do not use the equipment unless it is correctly earthed.

If you replace the plug keep in mind the following:


|                    |         |
|--------------------|---------|
| Blue wire:         | Neutral |
| Brown wire:        | Phase   |
| Yellow/green wire: | Ground  |

## 8 Operation

### STARTING UP

- Connect the equipment to the power supply.
- Turn on the main switch (12). The displays (1) and (2) will light.
- Press key (8) to open the lid.
- Insert the tube in the rotor in balanced compensate way.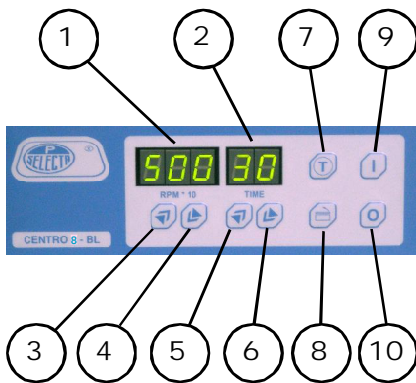

### HOW TO PLACE THE HEAD IN THE CENTRIFUGE CENTROLIT II-BL Placing the head code 7001704:


- Place the head in the axle and turn it until it fits perfectly with the bolt.
- Place the binding nut with the hexagonal side down and press as shown in photograph 1 and 2..
- Place the cover as shown in photo 3.

Placing the head code 7001703:

- Place the head in the axle similar as the previous head. See photo 4.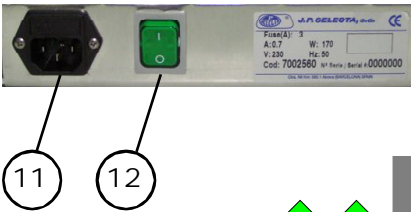

- Close the lid pressing on it until you heard the sound of the lock.
- By pressing the keys (3) and (4), you can increase or decrease the speed set which is shown on the display (1).
- By pressing the keys (5) and (6), you can increase or decrease the set time which is shown on the display (2).

**WORKING FOR A NOT DEFINED TIME**

When the display (2) shows «00» by pressing the key (6) the display shows: «--». If the centrifuge starts with this set time it will only stop manually.

- Start pressing the key (9). The centrifuge will stop once finished the set time.  
During the operation, the display (1) shows the speed read by the tachometer.  
The work can be stop by pressing the key (10).

**INSTANT START/STOP**

- Do the previous operation as explained on the section «Starting up». Once the speed of the operation is set and is different of 0, the rotor starts to turn at the speed set by pressing the key (7). When stop pressing, the centrifuge will stop.

**EMERGENCY OPENING**

In case of a cut off the power supply and if it would be necessary to open the door, please do the following:

- Remove the plastic stopper on the right lateral of the equipment.
- Pull carefully from this stopper and the lid will open.

**Do not use this emergency opening device when the machine is working.**

**ERRORS LIST:**

- Error 1: Eprom failure
- Error 3: Tachometer speed failure
- Error 4: Overcurrent
- Error 5: Low voltage
- Error 6: Motor feedback voltage too high

## 9 Maintenance

Before removing the cover to handle inside, unplug the equipment from the socket.

**The handle of the electronic circuit boards by unauthorized personnel may cause damages. Make sure to give the equipment to an authorised technical assistance.**

### 9.1 CLEANING

We recommend to use a cotton cloth wet with a water-alcohol disolution.

### 9.2 SPARE PARTS

The more common spare parts:

|                                 |         |
|---------------------------------|---------|
| Spare description               | 7002357 |
| Electromagnetic lock            | 10040   |
| Main switch                     | 20017   |
| Motor                           | 24226   |
| Tachometer sensor circuit board | 29230   |
| Control panel circuit board     | 29364   |
| Power circuit board             | 29367   |
| Mains filter                    | 45011   |
| Transformar 115-115/18V 1.2A    | 45048   |


## 10 Guarantee

This product is guaranteed for one year. The guarantee does not cover damages caused by incorrect use or causes beyond the control of J.P. SELECTA,S.A.

Any manipulation of the equipment by personnel not authorized by J.P. SELECTA,S.A. automatically cancels the guarantee.

# 11 Electric diagram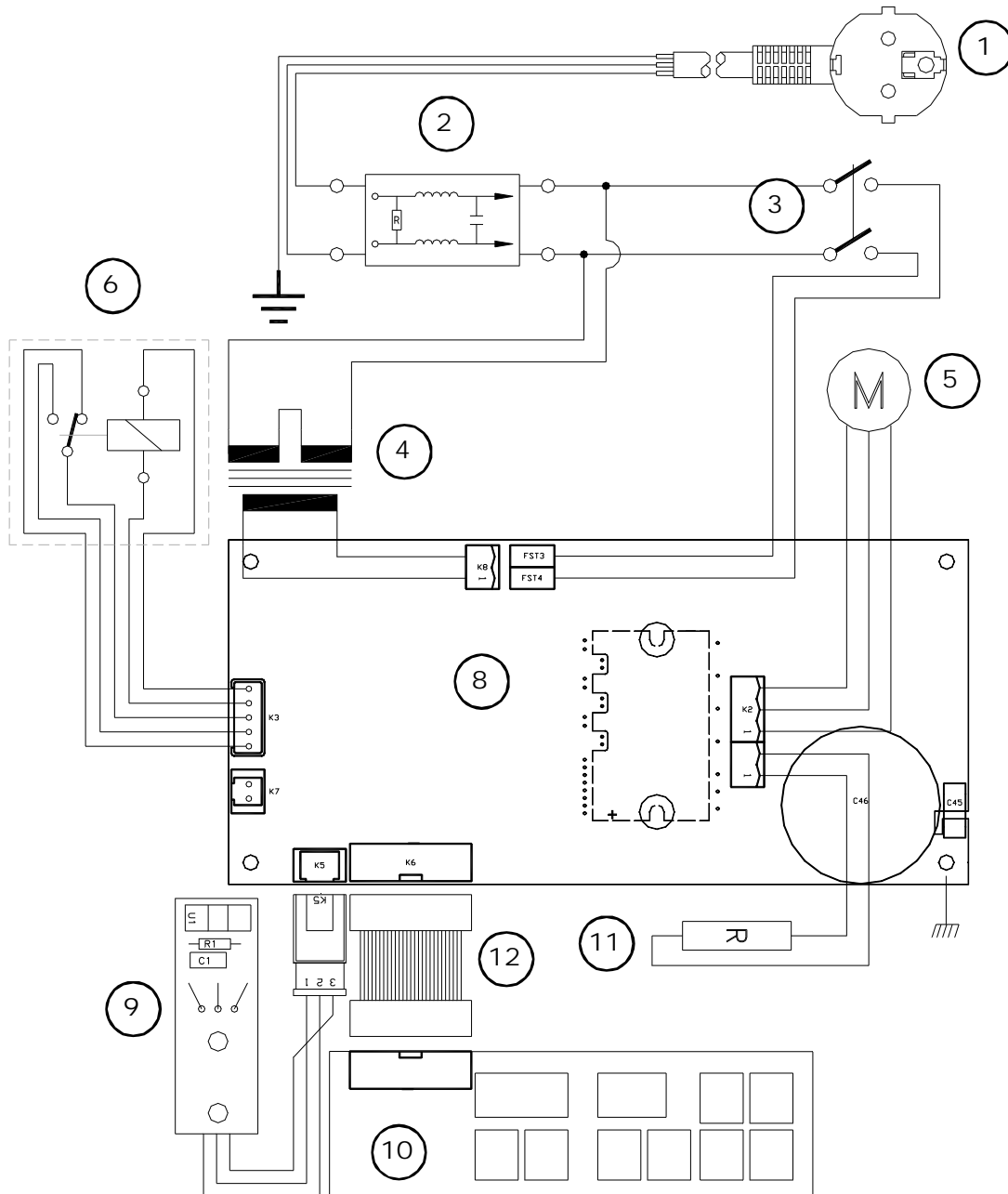

| Pos | Description                   | Code  |
|-----|-------------------------------|-------|
| 1   | Power supply cord             | 7140  |
| 2   | Mains filter                  | 45011 |
| 3   | Main switch                   | 29017 |
| 4   | Transformer 115-115/18V 1.2 A | 45048 |
| 5   | Motor                         | 24226 |
| 6   | Electromagnetic lock          | 10040 |

| Pos | Description                     | Code  |
|-----|---------------------------------|-------|
| 8   | Power circuit board             | 29367 |
| 9   | Tachometer sensor circuit board | 29230 |
| 10  | Control panel circuit board     | 29364 |
| 11  | Break resistance                |       |
| 12  | Board connecting wire           |       |

© J.P. SELECTA, s.a. 2007

---

---