CÓDIGO / PART NUMBER

1001403

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PRODUCT DESICCATOR

Indice

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English Version

1. General Information

- Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.
- Do not install or use the equipment without reading this handbook first. This instruction manual must be attached to the equipment and must be available for all users.
- If you have any doubts or enquiries, contact with J.P. SELECTA's technical service.

2. Packing List

Model:	1001403
Desiccator	1
Trays	3
Vessel for desiccant product (inox AISI 310)	1
Instructions manual	80273

3. Technical features

Capacity:	55
Inner dimensions: height/width/depth (cm)	50/38/29
Hygrometer (%HR)	0-100

4. Equipment description

Desiccator cabinet to store anhydrous, biologic or chemical products or samples which have to be preserved against humidity or duster.

It is made from transparent methacrylate. Door with a rubber silicone gasket and magnetic lock and a hydrometer graduated from 0-100% HR.

5. Installation and starting up

5.1 Unpacking

Once unpacked, visually check the equipment to see if it is in good condition. Keep the packaging box for some days.

5.2 Placing and assembling

1. Place the device on an even surface.
2. Assemble the three supplied shelves inside the cabinet.
3. Place the stainless steel tray on the bottom of the cabinet.

5.3 Desiccant product

Fill the desiccant product in the inner vessel. The recommended product is Silica Gel (SiO₂). It can be used with or without indicator. The desiccator with indicator allows to follow the degree of water absorption at first sight. The indicator of cobalt chloride looks blue when it is dry and changes to pink when it gets moisten.

6. Cleaning

The equipment can be cleaned with an alcohol or glass cleaner wet duster.

7. Warranty

This product has a warranty period of a year. The warranty does not include damages caused by an improper use or for any reason alien to J. P. SELECTA, s.a.u.

Any operation made by not authorized personnel cancels the warranty.

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