TEMPLOW-M

2101279

**ARMARIO CONGELADOR
UPRIGHT FREEZER**

Indice

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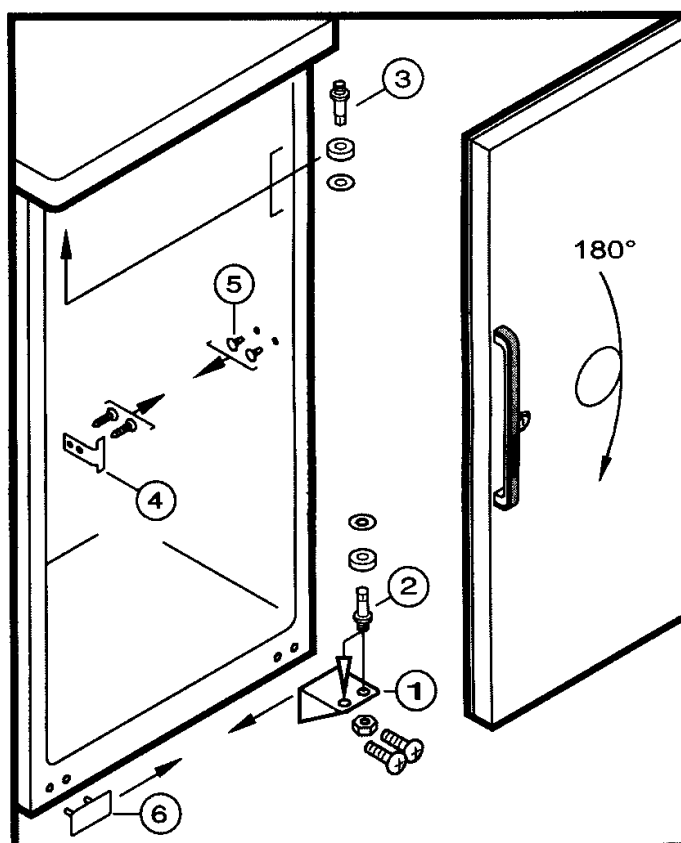
English version

1. GENERAL INFORMATION

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please inform the supplier immediately. To prevent injury or damage to the unit, it must be unpacked and set by two people.
- 2) If you get an equipment damaged on delivery, please contact the supplier immediately before connecting to the power supply.
- 3) Do not install or use the equipment without reading this handbook before.
- 4) To guarantee a safety operation, make sure that the equipment is set up and connected as described in this operating instructions.
- 5) When disconnecting the appliance, pull out from the plug, not from the cord.
- 6) This handbook must always be attached to the equipment and it must be available for all users.
- 7) Any doubt you have can be clarify by your supplier or the technical service of J.P. Selecta, s.a.u.
- 8) Any repairing or work on the appliance must only be carried out by the customer service.
- 9) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 10) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.
- 11) **DO NOT USE THE EQUIPMENT WITH FLUIDS WHICH CAN GIVE OFF VAPOURS CAPABLE OF MAKING EXPLOSIVE MIXTURES.**
- 12) **DO NOT ALLOW NAKED FLAMES OR IGNITON SOURCES INSIDE THE APPLIANCE.**
- 13) **DO NOT USE ELECTRICAL EQUIPMENT INSIDE THE APPLIANCE.**
- 14) **USE THE APPLIANCE IN A VENTILATED ROOM BECAUSE IN CASE OF REFRIGERANT LEAKAGE IT MIGHT CAUSE AN EXPLOSION.**

2. EQUIPMENT DESCRIPTION

The upright freezer **Templow-M** is useful to preserve products that require being stored at low temperatures. It can reach temperatures from -9°C until -26°C although the minimum temperature reached depends on the room temperature.

It is built of an epoxy covered external case and an internal ABS molded chamber which is high chemical resistant. And it also has rilsan coated steel baskets.

3. TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Capacity (litres):	520L		
Power (W):	345W		
Inner dimensions (cm):	Height 150	Width 64	Depth 53
External dimensions (cm):	Height 172	Width 75	Depth 72

4. INSTALLATION

- Avoid placing the appliance in direct sunlight or near cookers, heat radiators or similar heat sources.
- The floor on which the appliance stands should be horizontal and levelled. To compensate for any unevenness, please adjust the height of the legs.
- The ventilation grilles should not be obstructed. Always make sure that there is good ventilation and that the outward flowing air is able to escape.
- Do not place heat-emitting appliances on top of the freezer.
- Standard EN 378 specifies that the room in which you install your appliance must have a volume of 1m³ per 8g of R600a refrigerant used in the appliance, so as to avoid the formation of flammable gas/air mixtures in the room where the appliance is located in case of leakage in the refrigerant circuit. The quantity of the refrigerant used in your appliance is indicated on the type plate inside the equipment.

THE MINIMAL VOLUME IN THIS CASE IS 5 m³.

5. CONNECTION TO THE MAINS

Power supply and voltage at the operating point must meet the details on the features plate, which is located in the freezer compartment on the left side. The socket must be fused with a 10A fuse or higher, it must be away from the rear part of the appliance and must be easily accessible.

Connect the equipment with a properly earthed plug and socket.

If you have to change the plug cap, keep in mind the following:

Blue wire: Neutral.
Brown wire: Phase.
Yellow/green wire: Earth.

6. CONTROL PANEL

1. Temperature display.
2. Temperature buttons.
3. On/Off button.
4. Audible warning button.
5. Child-proof lock.

7. SWITCHING THE EQUIPMENT ON AND OFF

On/Off

Switching on: It is recommended to clean the appliance before switching it on for the first time (see "Cleaning" section). Press the **ON/OFF** button so that the temperature display lights up/flashes. When the appliance is switch on and while it is still warm, the display will contain dashes until the temperature falls below 0°C. Switch the appliance on 4 hours before loading it with frozen materials for first time.

Switching off: Press the **ON/OFF** button for approximately 2 seconds. The temperature display will switch off.

8. SETTING THE TEMPERATURE

Push **UP** button to increase temperature (less temperature) and **DOWN** button to reduce the temperature (more temperature).

- While you select the temperature, the value will be flashing.
- When you push the setting buttons for the first time, the most recent setting will be displayed.
- The setting value will be increased in increments of 1°C.
- After 5 seconds after having pressed the last button, real temperature will be displayed in the equipment.
- Temperature can be set between -9°C to -26°C.

9. TEMPERATURE DISPLAY

The thermometer shows the highest temperature of the frozen material. It is adjusted and designed so that only temperature below 0°C is displayed.

The temperature can rise temporarily because inflowing warm air when room temperature material is stored or when frozen goods are removed or rearranged.

If F3 appears in the display, the appliance has a system fault. Please contact the technical service and indicate this fault number.

10. WARNING SIGNAL

This equipment comes equipped with an alarm function.

- The warning signal sounds when the door has been left open for more than 60 seconds. Press the **ALARM** button to cancel it. When the door is shut, the warning signal is again in standby mode.
- When the freezer temperature is not cold enough, the warning signal will sound and the temperature display will flash. Press the **ALARM** button to cancel it. The temperature indication will continue flashing until the alarm situation has been rectified.

11. ADDITIONAL FUNCTIONS

There are two additional functions: adjusting brightness of display and child-proof lock, to avoid the machine being disconnected or modified accidentally.

Activate set-up mode:

- Press the **ALARM** button for about 5 seconds. The display will show a flashing **c**.
- Select the required function by pressing the **UP/DOWN** button.
 - c** = child-proof lock,
 - h** = display brightness.
- Press the **ALARM** button briefly to select/confirm.

Press  for 5 seconds. Display = 

 Display = 

 Display = 

 Display =  

 Display =  

If you select **c = child-proof lock**, push **UP/DOWN** button to change from **c0** (child-proof lock off) to **c1** (child-proof lock on) and confirm by pushing the **ALARM** button.

When the symbol  is lit, the child-proof lock is activated.

Press  for 5 seconds. Display = 

 Display = 

 Display = 

If you select **h = brightness**, push **UP/DOWN** button to change from **h1** (minimum) to **h5** (maximum brightness) and confirm by pushing the **ALARM** button.

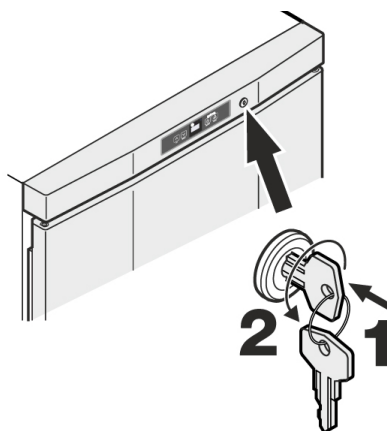
On/Off

Exit set-up mode:

To exit the set-up mode, push the **ON/OFF** button. After 2 seconds, the system will return to its normal operating mode.

12. SAFETY LOCK

The appliance lock is equipped with a safety mechanism. To lock the freezer, please insert the key in the 1 direction and turn it 90° to the left (2). To unlock the equipment, proceed in the same order.

13. DEFROSTING

After a long period of operation, a layer of frost or ice builds up on the cooling plates of the freezer compartment. This increases the energy consumption of the freezer. The cooling plates should therefore be defrosted regularly.

- To defrost, switch the appliance off. Pull out the mains plug.
- Wrap the frozen goods in paper (newspaper) or blankets and store them in a cool place (in the refrigerator, for instance).
- To speed up the defrosting process put a saucepan of hot (but not boiling) water on one of the cooling elements.
- Leave the door of the appliance open while defrosting. While defrosting, mop out the remaining water and clean the appliance.
- Do not use mechanic tools or other artificial aids to defrost as it may produce serious damages.

14. CLEANING

Before cleaning, disconnect the appliance from mains. Clean the equipment with warm water and a little detergent. Do not use abrasive or acid cleaners or chemical solvents, as they could damage the surface and cause corrosion. We recommend the use of a neutral pH cleaner. Do not use steam cleaners because of the risk of injury and damage.

- Empty the refrigerator and store the goods as it is indicated on the "Defrosting" section.
- Remove the storage shelves and support rails.
- Make sure that no cleaning water penetrates into the electrical components or ventilation grille.
- Do not damage or remove the type plate from the inside of the appliance. It is important for servicing purposes.
- Dry all parts well with a cloth, when finishing.
- Once a year, the dust should be removed from the refrigeration unit and the heat exchanger - metal grid at the back side of the appliance.

15. LONG TIME DISCONNECTION

If you must shut down the appliance for a long time period, switch it off and disconnect the plug. Also, clean it and leave the door open in order to prevent bad smells.

16. MALFUNCTIONS

In case of malfunction, you may be able to check the following:

- * **The appliance does not function, the control lamps do not lit:**
 - Is the appliance switched on?
 - Is the plug correctly fitted in the mains socket?
 - Is the fuse burned out?
- * **The temperature is not low enough:**
 - Is the temperature setting correct?
 - Have excessive amounts of material been placed in the appliance?
Check the displayed temperature after 24 hours.
 - Is the ventilation system working properly?
 - Is the appliance set up too close to a heat source?

If none of the above causes apply and you cannot rectify the fault yourself, please contact the technical service.

17. CHANGING OVER DOOR HINGES

- Close the door and unscrew the lower hinge bracket (1).
- Lift the door, tilt it to the right and remove it.
- Rotate 180° and remove the lower hinge spin (2) and screw it on the opposite side.
- Unscrew upper hinge elements (3) and screw them on the opposite side.
- Attach the hinge (4) and the plugs (5) to the opposite side.
- Transfer lower hinge bracket and cover plate (6) to the opposite side.
- Locate door on upper pin.
- Tilt door inwards from the left and secure the hinge pin with a screw.
- Transfer handle and plugs to the opposite side of the door.